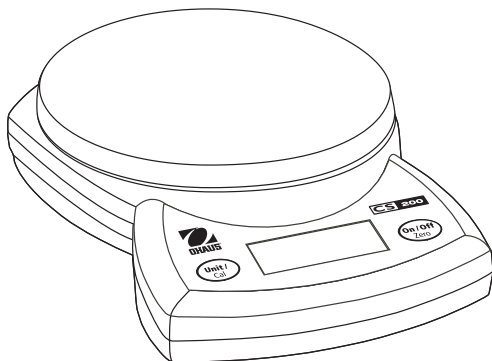
Ohaus Corporation
19A Chapin Road
P.O. Box 2033
Pine Brook, NJ 07058-2033, USA

Ohaus Compact Scales Models CS 200, CS 2000 and CS 5000 -OPERATING MANUAL-

- The CS 200, CS 2000 and CS 5000 only differ in weighing capacity and readability. Unless otherwise specified, the content of this manual applies to all three models.
- For best results, use the scale in the specified optimum operating temperature.
- For best accuracy, perform your mass measurements in clean environments, free of electrostatic and magnetic fields, dust and dirt, moisture, vibration and air currents.
- Use your scale with care, gently loading items to be weighed onto the center of the pan. Avoiding rough treatment will aid in the life of your compact scale.
- Calibration weights are not provided with the scale.

Battery Installation

Three AA size alkaline batteries are provided.

- Remove the battery cover on the bottom of the scale and place the three “AA” size alkaline batteries into the compartment as indicated.
- Do not use excessive force or press the weighing pan.
- Re-insert the battery cover.

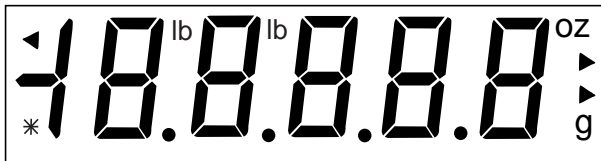
AC Adapter Operation

AC power adapters are available as optional accessories.

- Connect the AC adapter to the power input jack at the rear of the scale.
- Plug the AC adapter into a properly grounded wall outlet.

General Product Features

- **Power Up Test:** When the scale is turned on, all display segments will appear for approximately 3 seconds before resetting to zero.

- **Stable Reading Indication:** On the display, a star indicator will appear in the lower left corner of the display when a stable reading has been reached.
- **Weight Unit Indication:** On the display, either lb oz, oz, or g will appear indicating the weight unit selected.
- **Overload:** If the applied load exceeds the capacity of the scale, an E will appear on the display and the load should be removed immediately. The scale will return to normal operation.
- **Zero Function:** Values can progressively be added to a sample. By pressing the **On/Off Zero** button, the scale display returns to zero. After adding additional mass, press **On/Off Zero** to zero the scale again. Additional mass may be added up to the capacity of the scale.

- **Negative Value:** When a load is removed from the scale, any zeroed value will be displayed as a negative number. To return to normal operation, the zeroed value can be cancelled by pressing **On/Off Zero**.
- **Low Battery Indication:** The display will show Lo when the batteries are weak and need to be replaced.
- **Auto Shut-Off:** To extend battery life, the scale will automatically turn off after approximately five minutes if no button press or active weighing is occurring. This feature is only active with battery operation only.

Two Button Keypad

- **On/Off Zero:** Pressing this button turns on the scale. This same button operates the zero feature when the scale is on. Press and hold this button for three seconds to turn the scale off.
- **Unit/Cal:** Press this button briefly to change the weighing unit. The following weight units may be selected: pound-ounces (lb oz), ounces (oz), or grams (g). Press and hold this button to begin the calibration process.

Calibration

For best results, calibrate the scale at regular intervals. This is especially important if the scale is in use for prolonged periods.

- Pressing and holding the **Unit/Cal** button starts the calibration process. The display should show CAL.
- The calibration process can be aborted by turning the scale off.
- Press the **On/Off Zero** button to capture 0. The display shows -C- while the scale stores the the zero load signal.
- The display will show C xxx where xxx is the calibration weight in grams.
- Place the appropriate calibration weight on the platform.
- Press **On/Off Zero** button.
- The display shows -C- while the scale stores the calibration point signal.

A rectangular digital display showing the text "CAL" in a seven-segment font.

A rectangular digital display showing the text "-C-" in a seven-segment font.

A rectangular digital display showing the text "C 1000" in a seven-segment font.

A rectangular digital display showing the text "-C-" in a seven-segment font.

- After span capture, the display returns to the normal weighing mode.

- The message CAL E will appear if the calibration steps are not followed or the wrong weight was used.

Accessories Supplied

- Three AA size alkaline batteries.

Optional Accessories

AC Power Adapters:	Ohaus Part No.
120V US Plug	90524-66
220V Euro Plug	90524-63
240V UK Plug	90524-64
240V Australian Plug	90524-65
Calibration Masses:	
200g for CS 200	51025-16
1000g for CS 2000	51016-16
2000g for CS 5000	51026-16
1050 ml Bowl	80850075
Stainless steel weighing pan cover	80850076

Specifications

	CS 200	CS 2000	CS 5000
Capacity x Readability	200 X 0.1g 0 lbs: 7.05 oz x 0.01 oz	2000 X 1g 4 lbs: 6.55 oz x 0.05 oz	5000 x 2g 11 lbs: 0.4 oz x 0.1 oz
Linearity	±0.1g	±1g	±2g
Weighing units	g, lb, oz		
Calibration	User calibration		
Stabilization time	3 seconds		
Maximum overload	150% of capacity		
Overload indication	[E]		
Underload indication	Negative value		
Auto shut-off	5 minutes no activity		
Optimum operating temperature	64° to 77° F / 18° to 25° C		
Weigh pan size L x W (in/cm)	5.75 x 5.25/14.5 x 13.3		
Power requirements	3-AA alkaline batteries		
Dimensions L x W x H (in/cm)	8.00 x 5.50 x 1.50 / 20.5 x 14.0 x 3.8		
Net Weight (lb/kg)	1.0/0.47		


Declaration of Conformity We, Ohaus Corporation, declare under our sole responsibility that the balance models listed below marked with "CE" are in conformity with the directives and standards mentioned.

Scale model (s) Compact CS 200, CS 2000 and CS 5000

Marked with:	EC Directive (Including applicable amendments)	Standard
	73/23/EC Electrical equipment for use within specified voltage limits	EN60950:1992 Safety of Information Technology Equipment
	89/336/EC Electromagnetic compatibility	EN61326:1997 + A1:1998 Electrical Equipment for Measurement, Control and Laboratory Use - EMC requirements Immunity - Performance Criterion B Emission - Class B

Last two digits of the year which the CE marking was affixed: 02

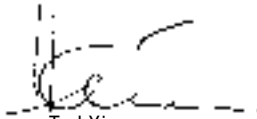ISO 9001 Registration for Ohaus Corporation. Ohaus Corporation, USA, was examined and evaluated in 1994 by the Bureau Veritas Quality International (BVQI) and was awarded ISO 9001 registration. This certifies that Ohaus Corporation, USA, has a quality system that conforms with the international standards for quality management and quality assurance (ISO 9000 series). Repeat audits are carried out by BVQI at intervals to check that the quality system is operated in the proper manner.

Ted Xia
President

Ohaus Corporation, Pine Brook, NJ USA

NOTE: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

LIMITED WARRANTY

Ohaus compact scales are warranted against defects in materials and workmanship for a period of 12 months from the date of delivery. During the warranty period, Ohaus will repair, or, at its option, replace any component(s) that prove to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

Ohaus Corporation
19A Chapin Road
P.O. Box 2033
Pine Brook, NJ 07058-2033, USA
Tel: (973) 377-9000
Fax: (973) 593-0359

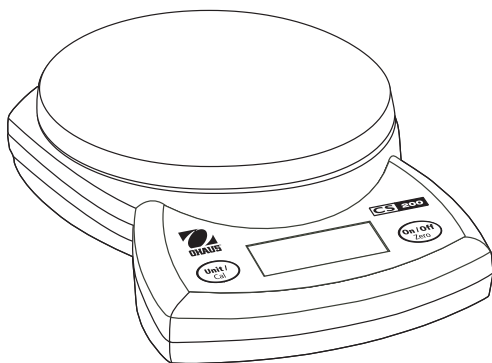
With offices worldwide.

Ohaus Corporation 2002, all rights reserved.

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