

Tray capacity & flask positioning


The TF18 flask tray contains threaded holes to hold 25ml, 50ml, 100ml, 250ml, 500ml and 1000ml flask clamps.

The following table and diagrams show the tray capacity and optimal clamp positions.


TF18	25ml	50ml	100ml	250ml	500ml	1000ml
Flask capacity	33	26	17	14	6	4

Attaching the clamps to the flask tray


- M4 fixings are supplied with all flask clamps:
 - 25ml, 50ml & 100ml clamps have one fixing
 - 250ml, 500ml and 1000ml clamps have four fixings
- Secure the flask clamps onto the tray using a suitable screwdriver

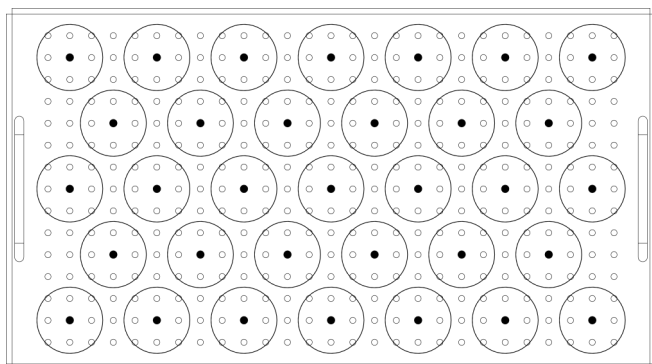


25ml

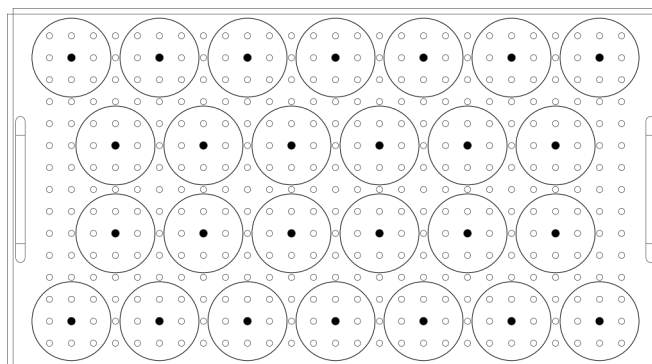


50ml

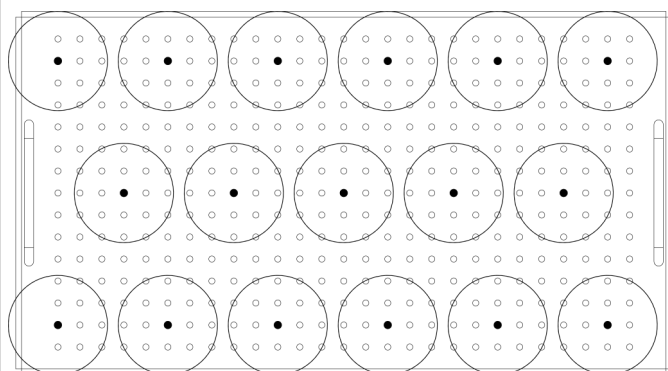


100ml

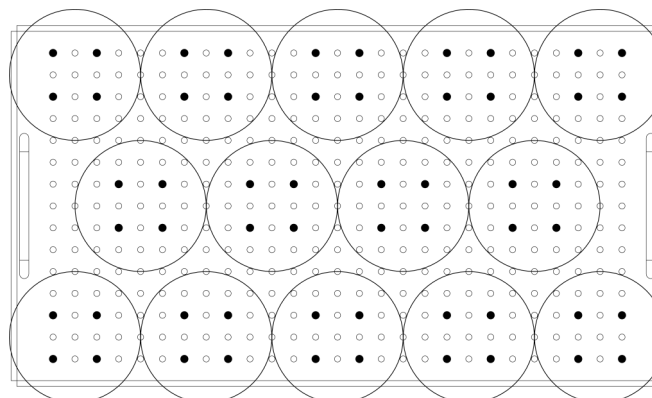

250ml

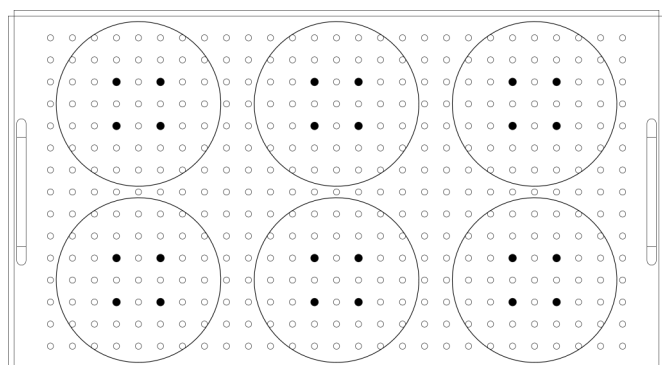

500ml

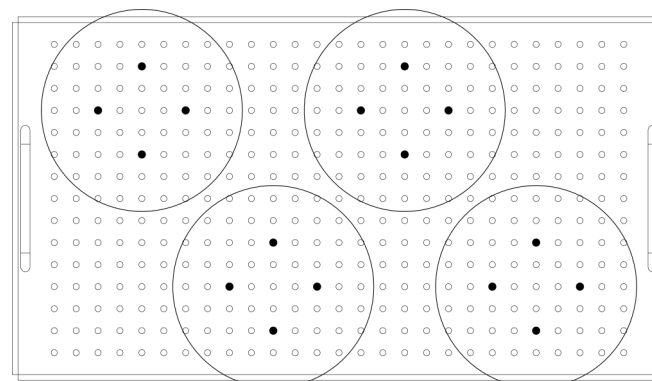

1000ml