Lovibond® Colour Measurement

Tintometer® Group





- √ Hand-Held Imaging Spectrocolorimeter
- √ Unique housing stand to protect sample from ambient light
- √ Enables measurement of food, powders, cosmetics & pastes
- √ Versatile Communication of Colour Measurement
- √ Single, Simple, Reliable, Integrated Tool no need to connect to PC
- √ Readings and Comparison Charts within Seconds
- √ Information on Colour and Surface Texture Effects
- √ Conformance Reports in PDF and Excel Formats
- √ Precise Image Sampling with 9 Band Accuracy

Lovibond® LC 100

The Lovibond® LC 100 enables setting and maintaining colour standards throughout the manufacturing process. It is a lightweight tool that is easy to handle and ready to use. In just over a second, it not only provides a reading of the sample but also a report telling how close the colour is to a measured standard. Together with the Lovibond® SV 100, it is the ideal solution for the colour measurement of food, powder, cosmetics, pastes and pharmaceuticals.

The LC 100 automatically creates conformance reports in either PDF or Excel formats for each stored sample against each standard. These can easily be shared with a PC or directly to a compatible printer (via USB mass storage mode).

Features

- Enhanced Working Agreement:
 Takes images of sample using 8 different visible illuminations and
 1 UV LED (9 bands) for more accurate colour measurement
- Unique Imaging Technology: 45/0 optical geometry and image capture technology: better agreement to visual observations
- Information on Colour and Surface Texture Effects:
 Each measurement is a combination of 27 images illuminated with different colours and from different directions
- Quick, Accurate Reading:
 - Measurement time of just over a second
- Precise Positioning on sample:
 - Enables preview of measurement area and storage and recall of image after measurement
- Increased Quality Control:
 - Up to 20 reference standards can be measured and stored each date and time stamped with option to name $\,$
- Single Unit for Multiple Tests:
 - Up to 350 sample measurements can be stored on the device and compared to the reference standards
- Easily Identify Measurements:
 - Samples can be tagged via the on-screen keyboard for text or via the built in voice recorder
- Save and Print Results:
 - For a more in-depth report, the LC 100 creates PDF and CSV files which can be accessed via a PC with the USB mass storage mode

Test Results

Test results are easily viewed with:

- √ Graphical Pass/Fail indicator
- \checkmark Display of ∆E for all common methods ie CIELAB, CMC, CIE 94, 2000
- √ Display of measured and delta differences with L*a*b*C*h° for standard and sample
- √ Graphical L*a*b* plot
- √ Verbal colour difference description



The Tintometer Ltd T: +44 (0)1980 664800 F: +44 (0)1980 625412 Tintometer North America T: +1 941 758 8671 F: +1 941 727 9654

Toll Free: 855-Color12 (855-265-6712)

Lovibond® SV 100

The Lovibond® LC 100 and SV 100 integrated package provides a unique solution to users looking to measure food, powders, cosmetics and pastes.

The SV 100 holder supports the LC 100 at an ergonomic viewing angle to enable the measurements in Lovibond® optical cells. A secure cap is provided to cover the sample, protecting it from any excessive ambient light, enabling quick yet accurate readings.

The units are provided in a purpose-designed case containing all the accessories required for accurate colour measurement expected of the Lovibond® brand.





Technical Specification

Measuring Geometrics 45/0 Image Capture

Light Source Independent tri-directional 25 LED (8* visible wavelengths; 1* UV)

Illuminant/Observer D65&A(10)

Standards / Sample Storage 20/350

Measurement Time 1.8 seconds

Measuring Area 4 & 8mm

Short Term Repeatability Typical 0.10 ΔE 94 on white

(D65/10)

Display 4.5cm Colour TFT

Data Interface USB (Mass Storage Device)

Operating Temperature 0° to 40°C (50° to 104°F)

Storage Temperature -20° to 60°C (-4° to 140°F)

Humidity Range 20-80% RH (non-condensing)

Usage Indoor Use Only

Altitude 2000m

Polution Degree 2

Transient Overvoltage Category II

	Order Number	Product
	401100	Lovibond® LC 100 Spectrophotometer
	401120	Integrated Package : LC 100 & SV 100 Kit
	132750	Replacement Cells (10mm, Pk 100)

Tintometer - South East Asia T: +603 3325 2286 F: +603 3325 2287 Tintometer - China T: +86 10 8525 1111 Ext 330 F: +86 10 85251001

sales@tintometer.com