

REF 918 073

04.12

en

NANOCOLOR[®] Chloride detection kit AOX

Method:

Precipitation test with silver nitrate

Contents:

10 mL silver nitrate solution 1–5 %

Hazard warning:

For further information please ask for a safety data sheet.

Procedure:

After rinsing, as described in test 0-07 AOX 3, proceed as follows:

Rinse NANOSORB cartridge with
5 mL distilled water and collect the water in a clean beaker or tube.

Add

3 drops of silver nitrate solution and mix.

Solution clear: go on with test 0-07 AOX 3

Solution turbid: rinse again with rinsing solution

Disposal:

The contents of the beaker can be washed into drain with plenty of water.

MACHEREY-NAGEL GmbH & Co. KG

Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

Tel.: +49 24 21 969-0 · Fax: +49 24 21 969-199 · info@mn-net.com · www.mn-net.com

918073/04.2012