

Incubating Cooling Orbital Shakers



OHAUS Incubating Cooling Orbital Shakers are designed to incubate samples from 10°C below ambient to 65°C. Incubating Cooling Mini Shaker is designed to hold two microplates or two optional modular tube blocks in a variety of configurations. All models feature touchpad controls with easy-to-read, independent LED displays for temperature, speed and time. Microprocessor control provides consistent uniform shaking action.

- Microprocessor-Controlled for Consistent Shaking Action
- Triple Eccentric Drive Provides Reliable Service and Continuous Duty
- Safety Features Include Speed Ramping and Overload Protection

Incubating Cooling Orbital Shakers

- Heats to 65°C and cools to 10°C below ambient
- LED displays for temperature, speed and time
- Calibration mode for temperature

The OHAUS Incubating/Cooling Orbital Shaker is microplate ready without the need for any additional accessories. Optional modular blocks can accommodate micro-tubes, centrifuge tubes, vials, or culture tubes. Unit holds microplates or modular blocks with a 12.7 cm tall interior capacity. Ideal for analyses that require a stable, controlled temperature.

Operating Features:

Microprocessor Control: The variable speed microprocessor control provides consistent uniform shaking action. Microprocessor will display last set-point and will restart if power is interrupted.

PID Temperature Controller: Maintains precise temperature control from 10°C below ambient to 65°C. Easy-to-use controls allow users to adjust temperature in 1°C increments.

Triple Eccentric Drive: Permanently lubricated ball bearings and maintenance free, brushless DC motor provide reliable service and continuous duty operation.

LED Display: Touch pad controls with easy-to-read, independent LED displays for temperature, speed, and time allow operator to view all settings at once. Provides repeatable and accurate results every time and is easily visible across lab benches. Timer will display elapsed time or, when programmed to user defined limit, will shut off unit when time reaches zero. Display will show last used settings, even after power has been turned off.

Temperature Calibration Mode: Allows user to calibrate unit to an external temperature device.

Safety Features:

Overload Protection: Audible and visual signals will activate when system detects an obstruction or overload of the tray.

Speed Ramping Feature: Slowly increases speed to desired set-point to avoid splashing.

Audible Alarm: In timed mode, alarm will sound when the time reaches zero. Alarm has optional mute function which can be set from the touch pad control.

Caution Hot Indicator: Symbol illuminates when the temperature of the air in the chamber reaches 40°C and remains lit until temperature cools down.

Spill-Resistant Design: Channels fluids away from internal components.

Polycarbonate Lid: Permits viewing of samples without disturbing internal temperature.

Operating Conditions:

Unit can be run in conditions from 5 to 40°C, maximum 80% relative humidity, non-condensing.

Applications:

Cell and bacterial cultures, hybridizations, and enzyme reactions.

Ordering Information:



Unit includes a detachable, 3-wire cord and plug. Units are also supplied with an adapter bracket to hold optional modular blocks. See page 36-37 for block options.



Specifications	
Temperature Range	10°C below ambient to 65°C
Temperature Uniformity	± 0.5°C at 37°C
Speed Range using Microplates	100 to 1200 rpm
Speed Range using Modular Blocks	100 to 600 rpm
Speed Accuracy	± 2%
Timer	1 second to 160 hours
Orbit	3 mm
Maximum Weight Capacity	2 microplates or 2 modular blocks
Overall Dimensions (L × W × H)	45.5 × 27.9 × 26.7 cm
Ship Weight	15.4 kg

Description	Model	Item Number
Incubating Cooling Light Duty Shaker	ISICMBCDG	30391940

Incubating Light Duty Orbital Shakers Modular Blocks

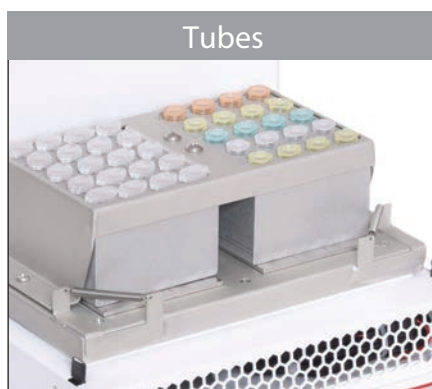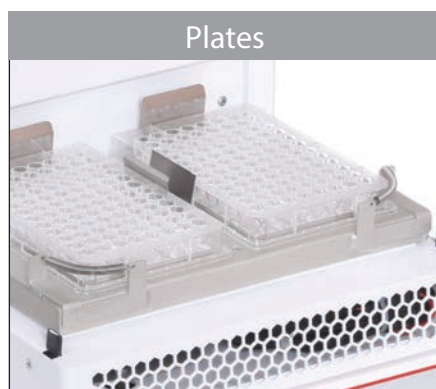

Modular Blocks

Modular Blocks are constructed from a solid anodized aluminum block. The close contact of tubes to block walls allow for maximum temperature transfer.

Block dimensions (L × W × H): 9.5 × 7.6 × 5.1 cm

Applications: Cell cultures, hybridizations, and extraction procedures


OHAUS modular blocks also fit in the OHAUS Dry Block Heaters.

Constructed of anodized aluminum, this material is ideal for its temperature conducting and corrosion resistant properties.

Microcentrifuge Tube Blocks

Single block

Sample Type	No. of Wells	Well Diameter	Well Depth	Item Number
0.5 mL tube	30	7.9 mm	27.6 mm	30400157
1.5 mL tube	20	11.1 mm	39.1 mm	30400159
2 mL tube	20	11.5 mm	38.1 mm	30400191

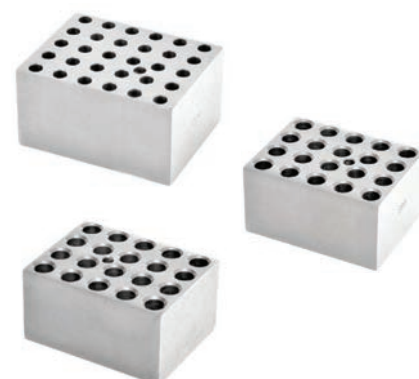

Conical-Bottom Centrifuge Tube Blocks

Single block

Sample Type	No. of Wells	Well Diameter	Well Depth	Item Number
15 mL tube	12	17.1 mm	44.5 mm	30400172
50 mL tube	5	29.0 mm	47.6 mm	30400168



Incubating Light Duty Orbital Shakers Modular Blocks

Standard Test Tube Blocks

Single block

Sample Type	No. of Wells	Well Diameter	Well Depth	Item Number
6 mm tube	30	8.3 mm	48.4 mm	30400158
10 mm tube	24	10.7 mm	48.4 mm	30400151
12/13 mm tube	20	13.9 mm	48.4 mm	30400152



Combination Blocks

Single block

These blocks have been designed for variable sized samples.

Sample Type	No. of Wells	Well Diameter	Well Depth	Item Number
Test Tube Combination				
6 mm	6	8.3 mm	48.4 mm	30400156
12/13 mm	5	13.8 mm	48.4 mm	
25 mm	3	26.2 mm	48.4 mm	
Centrifuge Tube Combination				
1.5 mL	4	11.1 mm	39.1 mm	30400193
15 mL	3	17.1 mm	44.5 mm	
50 mL	2	29.0 mm	47.6 mm	
Micro-Tube Combination				
0.5 mL	6	7.9 mm	27.6 mm	30400194
1.5 mL	10	11.1 mm	39.1 mm	
2 mL	5	11.5 mm	38.1 mm	



Vial Blocks

Single block

Designed for sample/serum and scintillation vials.

Sample Type	No. of Wells	Well Diameter	Well Depth	Item Number
12 mm vial	20	12.7 mm	30 mm	30400182
15 mm vial	20	15.8 mm	35 mm	30400183
16 mm vial	15	16.4 mm	45 mm	30400190
17 mm vial	12	17.8 mm	45 mm	30400184
19 mm vial	12	19.7 mm	45 mm	30400185
21 mm vial	9	21.7 mm	45 mm	30400186
23 mm vial	8	23.8 mm	45 mm	30400187
25 mm vial	8	25.8 mm	45 mm	30400188
28 mm vial	6	28.8 mm	45 mm	30400189

