

# *Incubating Light Duty Orbital Shakers*



OHAUS Incubating Light Duty Shakers are designed to incubate samples from 10° below ambient to 65°C depending on the model. The Incubating Mini Shaker has an 3.6 kg capacity while the Incubating Microplate Shaker can hold up to four standard or deep well plates. Incubating-Cooling Mini Shaker is designed to hold two microplates or two optional modular tube blocks. All models feature LED displays with touchpad controls.

- Microprocessor-Controlled for Consistent Shaking Action
- Triple Eccentric Drive Provides Reliable Service and Continuous Duty
- Safety Features Include Speed Ramping and Overload Protection

# Incubating & Incubating Cooling Shakers

## Incubating Light Duty Orbital Shakers

- LED displays for temperature, speed and time
- Timer with audible alarm
- Calibration mode for temperature

The OHAUS Incubating Light Duty Orbital Shakers are designed to heat and shake a variety of samples.

### Operating Features:

**Microprocessor Control:** The variable speed microprocessor control provides consistent uniform shaking action. Microprocessor will display last set-point and will restart if power is interrupted.

**PID Temperature Controller:** Maintains precise temperature control from ambient +5°C to 65°C. Easy-to-use controls allow users to adjust temperature in 1°C increments.

**Triple Eccentric Drive:** Permanently lubricated ball bearings and maintenance free, brushless DC motor provide reliable service and continuous duty operation.

**LED Display:** Touch pad controls with easy-to-read, independent LED displays for temperature, speed and time allow operator to view all settings at once. Provides repeatable and accurate results every time and is easily visible across lab benches. Timer will display elapsed time or, when programmed to user defined limit, will shut off unit when time reaches zero. Display will show last used settings, even after power has been turned off.

**Temperature Calibration Mode:** Allows user to calibrate unit to an external temperature device.

### Safety Features:

**Overload Protection:** Audible and visual signals will activate when system detects an obstruction or overload of the tray.

**Speed Ramping Feature:** Slowly increases speed to desired set-point to avoid splashing.

**Audible Alarm:** In timed mode, alarm will sound when the time reaches zero.

**Caution Hot Indicator:** Symbol illuminates when the temperature of the air in the chamber reaches 40°C and remains lit until temperature cools down.

**Spill-Resistant Design:** Channels fluids away from internal components.

**Polycarbonate Lid:** Permits viewing of samples without disturbing internal temperature.

### Operating Conditions:

Unit can be run in conditions from 5 to 40°C, maximum 80% relative humidity, non-condensing.

### Applications:

Cell cultures, bacterial suspensions, and hybridizations.

### Ordering Information:

Unit includes a detachable, 3-wire cord and plug.



Specifications	
Temperature Range	Ambient +5°C to 65°C
Temperature Uniformity	± 0.5°C at 37°C
Speed Range	100 to 1200 rpm
Speed Accuracy	± 2% of set speed
Timer	1 second to 160 hours
Orbit	3 mm
Maximum Weight Capacity	3.6 kg
Tray Material	Aluminum
Tray Dimensions (L × W)	27.9 × 19.7 cm
Interior Dimensions (L × W × H)	28.7 × 21.1 × 14.5 cm
Overall Dimensions (L × W × H)	43.2 × 27.9 × 27 cm
Ship Weight	13.6 kg

Description	Model	Item Number
Incubating Light Duty Orbital Shaker	ISLD04HDG	30391919

## Incubating Light Duty Orbital Shakers

- LED displays for temperature, speed and time
- Timer with audible alarm
- Available with opaque lid for light sensitive samples

The OHAUS Incubating Microplate Shakers are optimized for shaking microplates, deep-well plates, or micro-tubes.

### Operating Features:

**Microprocessor Control:** The variable speed microprocessor control provides consistent uniform shaking action. Microprocessor will display last set-point and will restart if power is interrupted.

**PID Temperature Controller:** Maintains precise temperature control from ambient +5°C to 65°C. Easy-to-use controls allow users to adjust temperature in 1°C increments.

**Triple Eccentric Drive:** Permanently lubricated ball bearings and maintenance free, brushless DC motor provide reliable service and continuous duty operation.

**LED Display:** Touch pad controls with easy-to-read, independent LED displays for temperature, speed, and time allow operator to view all settings at once. Provides repeatable and accurate results every time and is easily visible across lab benches. Timer will display elapsed time or, when programmed to user defined limit, will shut off unit when time reaches zero. Display will show last used settings, even after power has been turned off.

**Temperature Calibration Mode:** Allows user to calibrate unit to an external temperature device.

### Safety Features:

**Overload Protection:** Audible and visual signals will activate when system detects an obstruction or overload of the tray.

**Speed Ramping Feature:** Slowly increases speed to desired set-point to avoid splashing.

**Audible Alarm:** In timed mode, alarm will sound when the time reaches zero.

**Caution Hot Indicator:** Symbol illuminates when the temperature of the air in the chamber reaches 40°C and remains lit until temperature cools down.

**Spill-Resistant Design:** Channels fluids away from internal components.

**Polycarbonate Lid:** Clear lid permits viewing of samples without disturbing internal temperature. Opaque lid prohibits light exposure to light sensitive samples.

### Operating Conditions:

Unit can be run in conditions from 5 to 40°C, maximum 80% relative humidity, non-condensing.

### Applications:

Immunoassays and hybridizations.

### Ordering Information:

Units include a detachable, 3-wire cord and plug.



Specifications	
Temperature Range	Ambient +5°C to 65°C
Temperature Uniformity	± 0.5°C at 37°C
Speed Range	100 to 1200 rpm
Speed Accuracy	± 2%
Timer	1 second to 160 hours
Orbit	3 mm
Maximum Weight Capacity	4 microplates or 2 micro-tube racks
Tray Material	Aluminum
Tray Dimensions (L x W)	27.9 x 19.7 cm
Overall Dimensions (L x W x H)	43.2 x 27.9 x 19.7 cm
Ship Weight	13.6 kg

Description	Model	Item Number
Incubating Microplate Shaker	ISLDMPHDG	30391933
Incubating Microplate Shaker with Opaque Lid	ISLDMPHDGL	30391926