

Instructions for Mini EM4 Incubator

Supplied with:

- Handle – to be attached to the lid
- A perforated shelf
- Control thermometer on support

Specifications:

- Ideal room temperature for use: 20°C
- Temperature adjustable from ambient +5°C up to 44°C
- Useful volume: 4 litres
- External dimensions; 23cm W x 23cm D x 18cm H
- Internal dimensions; 18cm W x 18cm D x 12.5cm H
- Weight: 1.5kg
- Power: 25W, 220V

Operation:

- 1 - Temperature adjusting screw
- 2 - Power switch
- 3 – Red LED indicator – illuminates when the unit is hot



- Thermometer on support should be positioned in the centre of the incubator on the perforated shelf
- Do not put water into the incubator, doing so will void the warranty