

Customer Information

March 2020, LAW

Blanks in NANOCOLOR® Tube Tests

1. Required blank solutions for all NANOCOLOR® Photometers and therefore will be kept in the kits:

963026	COD Hg-free
985007	AOX 3
985015	Carbonate Hardness 15
985075	TOC 30
985078	TOC 300
985098	Aluminum 07

2. Water Zero. These tests can be measured on the PF-11 against the sample, water or the calibration cuvette:

985017	Chlorine/Ozone 2
985018	Chlorinedioxide 5
985024	Chromate 5
985031	Cyanide 08
985037	Iron 3
985043	Hardness 20
985054	Copper 7
985055	<i>ortho-</i> and <i>total</i> -Phosphate 45
985058	Manganese 10
985061	Nickel 7
985068	Nitrite 2
985069	Nitrite 4
985070	POC 200
985073	Sulfide 3
985074	Phenolic-Index 5
985076	<i>ortho-</i> and <i>total</i> -Phosphate 1
985080	<i>ortho-</i> and <i>total</i> -Phosphate 15
985081	<i>ortho-</i> and <i>total</i> -Phosphate 5
985084	Residual Hardness 1



3. Tests that need to be measured against a reagent blank:

985003	Ammonium 3
985004	Ammonium 10
985005	Ammonium 50
985006	Ammonium 200
985008	Ammonium 100
985012	COD 60 000
985019	Chloride 200
985021	Chloride 50
985023	COD 10 000
985026	COD 160
985028	COD 15 000
985029	COD 1500
985032	Detergents, anionic
985033	COD 300
985034	Detergents, cationic
985035	DEHA 1
985040	Fluoride 2
985041	Formaldehyde 8
985045	Potassium 50
985047	Detergents, nonionic
985050	Organic Acids 3000
985056	Molybdenum 40
985064	Nitrate 50
985079	<i>ortho</i> - and <i>total</i> -Phosphate 50
985083	<i>total</i> -Nitrogen TN _b 22
985085	Starch 100
985088	<i>total</i> -Nitrogen TN _b 220
985089	Sulfite 10
985091	Thiocyanate 50
985096	Zinc 4
985097	Tin 3

Contact

If you have further questions, please do not hesitate to contact us:

Technical support and customer service

Tel.: +49 24 21 969-332

E-Mail: csc@mn-net.com

For orders (international)

E-Mail: sales@mn-net.com

www.mn-net.com



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

DE/international:

Tel.: +49 24 21 969-0

Fax: +49 24 21 969-199

E-mail: info@mn-net.com

CH:

Tel.: +41 62 388 55 00

Fax: +41 62 388 55 05

E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68

Fax: +33 388 51 76 88

E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984

Fax: +1 484 821 1272

E-mail: sales-us@mn-net.com