

Customer Information

9th of June 2020, Version 3, RaKr

Firmware update for filter photometers NANOCOLOR[®] 500 D, 400 D und PF-12

Overview	
Photometer	Filter photometers NANOCOLOR [®] 500 D, 400 D und PF-12
Firmware version	2.71 (NANOCOLOR [®] 500 D) 3.70 (NANOCOLOR [®] 400 D) 2.69 (Compact photometer PF-12)
Price	Free of charge
Availability	From 09 th of June 2020
Available on	https://www.mn-net.com/nanocolor-device-updates
Required accessories	<ul style="list-style-type: none"> Update software for filter photometers Connection cable for the respective device



Important notes

The firmware update for the filter photometers NANOCOLOR[®] 500 D, 400 D and PF-12 erases the memory of the device. Make sure to save data prior to the update with our PC-software.

Update process

1. Unpack the file "Photometer-update-software-500D-400D-PF-12.zip" contained in this archive into a directory of your choice.
2. Open the file "NANOCOLOR@ Update.exe" to start the Firmware Update Software. You can change the language by choosing your desired language from the upper menu bar in the software.
3. Connect your photometer to your computer with the suitable connection cable for the respective device.
4. Turn on your photometer and type in method 900 to enter the settings.
5. Use the arrow keys of your photometer to navigate to "PROG. UPDATE". Confirm with the OK key.
6. In the sub menu pick "PROG. UPDATE" again and confirm with the OK key.
7. The firmware version ready to be installed on your photometer is shown in the Firmware Update Software on your computer.
8. Click on the button "Start update". The software automatically scans for the connected photometer and performs the firmware update.
9. After a completed update the message "Fotometer Update OK!" appears and you can now use your device with the most recent firmware.

Updating the update files

1. Click on „Internet" in the upper bar of the Firmware Update Software to automatically download the latest firmware version from our servers.
2. A popup informs, whether new firmware versions are available.

If this is not working, you can manually update the update files. Proceed as follows:

1. Unpack the file "Update-file-500D-400D-PF-12.zip" contained in this archive into the "data" folder of the Firmware Update Software directory.
2. Replace the files in the target directory.

www.mn-net.com



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

DE/international:

Tel.: +49 24 21 969-0
Fax: +49 24 21 969-199
E-mail: info@mn-net.com

CH:

Tel.: +41 62 388 55 00
Fax: +41 62 388 55 05
E-mail: sales-ch@mn-net.com

FR:

Tel.: +33 388 68 22 68
Fax: +33 388 51 76 88
E-mail: sales-fr@mn-net.com

US:

Tel.: +1 484 821 0984
Fax: +1 484 821 1272
E-mail: sales-us@mn-net.com

3. The most recent firmware versions are displayed in the software.
4. Perform the update as described in the chapter "Update process".

Changelog

- Implementation of new MN tests: *NANOCOLOR*[®] Manganese LR (REF 918126)
- Bugfixes

Contact

If you have further questions, please do not hesitate to contact us:

Technical support and customer service

Tel.: +49 24 21 969-332

E-Mail: csc@mn-net.com

For orders (international)

E-Mail: sales@mn-net.com

