



memmert
Experts in Thermostatics

Incubators

PERFECTLY COORDINATED. PERFECTLY CONTROLLED.



INCUBATOR I

CO₂ INCUBATOR IC_{Med}

COMPRESSOR-COOLED INCUBATOR IC_{Peco}/ICP

PELTIER-COOLED INCUBATOR IPP

COOLED STORAGE INCUBATOR IPS

100% ATMOSAFE. MADE IN GERMANY.

www.memmert.com | www.atmosafe.net



Stable. Safe. Sensitive.

Memmert incubators for microbiology.
Energy efficient, precise, 100% AtmoSAFE.

Even slight temperature deviations in the working chamber of an incubator may cause a test to fail. For this reason, the heating and control system of Memmert incubators are perfectly adapted to each other. During heating up and cooling down as well as in running operation, all appliances precisely keep the desired parameters within the smallest tolerance limits. Not only at one measuring point, but in the entire working chamber. Each individual Memmert incubator complies with the strict requirements of DIN 12880:2007-05 and is equipped with a maximum of safety functions. Each individual Memmert incubator is 100% AtmoSAFE.



INCUBATORS I

PAGE 4 - 8

Microbiological tests, colony counts, virology, toxicology

CO₂ INCUBATORS ICOmed

PAGE 9 - 13

Medical device class IIa for cultivation of cells or tissue, in-vitro fertilisation, gene expression

COMPRESSOR-COOLED INCUBATORS ICPeco

PAGE 14 - 18

Microbiological tests, colony counts, virology, toxicology, cultivation above and below room temperature, alternate stability tests

COMPRESSOR-COOLED INCUBATORS ICP

PAGE 19 - 20

Microbiological tests, colony counts, virology, toxicology, cultivation above and below room temperature, alternate stability tests

PELTIER-COOLED INCUBATORS IPP

PAGE 21 - 25

Protein crystallography, microbiological tests, colony counts, virology, toxicology, cultivation above and below room temperature, alternate stability tests

COOLED STORAGE INCUBATORS IPS

PAGE 26 - 29

Microbiological tests, cultivation above and below room temperature, stability tests

ADDITIONAL INFORMATION

PAGE 30





Incubator IN/INm and IF/IFm with SingleDISPLAY
Incubator INplus/INmplus and IFplus/IFmplus
with TwinDISPLAY
Natural convection or forced air circulation
AtmoCONTROL software

Model sizes:
30 / 55 / 75 / 110 / 160 / 260 / 450 / 750
+30 °C to +80 °C

INCUBATOR I Memmert incubators I are at home in the world of research, medicine, pharmaceuticals and food technology. Organic chamber loads require gentle heating. For this reason, the heating and control system are especially optimised for low temperatures of up to +80 °C. To prevent temperature overshoots, temperature is increased within a very narrow control range and kept exactly at the setpoint value. As required, the models with natural convection or with forced air circulation are available.



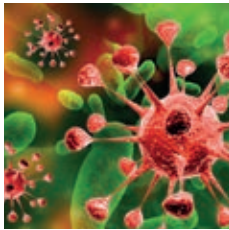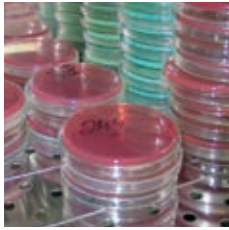

As little air circulation as possible in the incubator


Forced air convection may destroy the protective layer from moist air that is generated during incubation over the samples. This would lead to dehydration of the culture. In a Memmert incubator, the perfect combination of all-round surface heating and temperature control system ensures that incubation generally takes place without forced air circulation. Provided the chamber is fully loaded and forced air circulation is required, it can be precisely adjusted in 10 % steps from 0 to 100 %.

Sterilisation

The chamber of the incubators INplus/IFplus/INmplus and IFmplus, including all installations and sensors, can be sterilised at +160 °C in a 4-hour programme to guarantee optimum hygiene.

Fresh air is preheated

Temperature deviations caused by fresh air can influence sample characteristics or prolong drying. In Memmert incubators, the fresh air is therefore fed through a pre-heating chamber and seamlessly introduced into the working chamber.

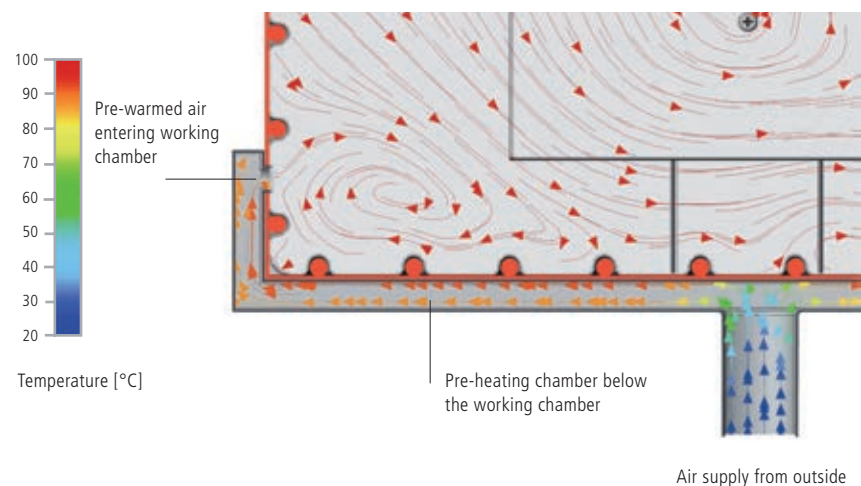

The incubator Im is a medical device:

The incubator Im is a Class I medical device in accordance with the EU directive 93/42/EEC. In accordance with the intended use incubators INmplus and IFmplus may be used for warming of rinsing solutions and infusions as well as of contrast agents. Incubators INm (with option A6) are intended for heating fango, silicate and APS packs for physical therapy and keeping them warm.

INCUBATORS I

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:
(EAC not valid for medical devices)



Interior: Stainless steel, material 1.4301 (ASTM 304) with all-round deep-drawn ribs to integrate the large-area heating with ceramic-metal sheath

Housing: Textured stainless steel, rear zinc-plated steel, intuitively operated SingleDISPLAY or TwinDISPLAY (TFT colour display) with touchscreen; inner glass door, outside fully insulated stainless steel door (from size 450 two leaves)


Fresh air: Admixture of pre-heated fresh air by electronically adjustable airflow

Connection: Mains cable with plug (German type)

Installation: 4 feet; sizes 450/750 mounted on lockable castors

Interfaces: Ethernet LAN, USB

USB: only TwinDISPLAY

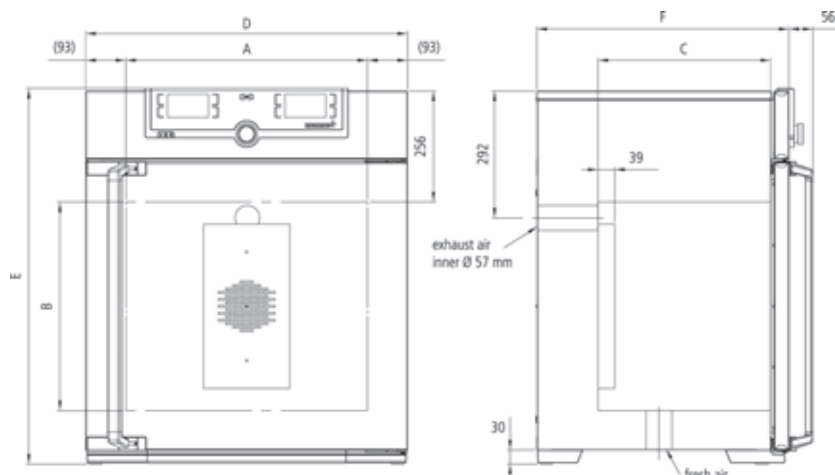

Model sizes/Description			30	55	75	110	160	260	450	750
Stainless steel interior	Volume	approx. l	32	53	74	108	161	256	449	749
	Width	(A) mm		400		560		640	1040	
	Height	(B) mm	320	400	560	480	720	800	720	1200
	Depth (less 39 mm for fan)	(C) mm	250	330		400		500	600	
	Max. number of grids/shelves	number	3	4	6	5	8	9	8	14
	Max. loading per grid/shelf	kg			20				30	
	Max. loading of chamber	kg	60	80	120	175	210		300	
	Max. loading per slide-in drip tray	kg		1,5		3		4	8	
	Max. loading per bottom drip tray	kg		1,5		3		4	8	
Textured stainless steel exterior	Width	(D) mm		585		745		824	1224	
	Height (size 450, 750 with castors)	(E) mm	704	784	944	864	1104	1183	1247	1726
	Depth (without door handle), door handle + 56 mm	(F) mm	434	514		584		684	784	
Standard equipment	Stainless steel grids, electropolished	number		1			2			
	Standard works calibration certificate (measuring point chamber center)	°C					+37			
Temperature	Working temperature range	°C	min. 5 (IN/INplus/INm/INmplus) 10 (IF/IFplus/IFm/IFmplus) above ambient temperature up to +80							
	Setting temperature range	°C	+20 to +80							
	Setting accuracy	°C	0.1							
Further data	Electrical load at 230 V, 50/60 Hz	approx. W	1600	1000	1250	1400	1600	1700	1800	2000
	Electrical load at 115 V, 50/60 Hz	approx. W	800			900			1500	1800
Packing data	Net weight	approx. kg	48	57	66	76	96	110	161	217
	Gross weight (packed in carton)	approx. kg	64	76	85	101	122	161	227	288
	Width	approx. mm	660	730		830		930	1330	
	Height	approx. mm	890	950	1130	1050	1300	1380	1440	1910
	Depth	approx. mm	650	670		800		930	1050	

Order No. Incubators

I = Incubator

N = Natural convection

F = Forced convection

m = Medical device

plus = Model with TwinDISPLAY

IN30 IN30m	IN55 IN55m	IN75 IN75m	IN110 IN110m	IN160 IN160m	IN260 IN260m	IN450 IN450m	IN750 IN750m
IN30plus IN30mplus	IN55plus IN55mplus	IN75plus IN75mplus	IN110plus IN110mplus	IN160plus IN160mplus	IN260plus IN260mplus	IN450plus IN450mplus	IN750plus IN750mplus
IF30 IF30m	IF55 IF55m	IF75 IF75m	IF110 IF110m	IF160 IF160m	IF260 IF260m	IF450 IF450m	IF750 IF750m
IF30plus IF30mplus	IF55plus IF55mplus	IF75plus IF75mplus	IF110plus IF110mplus	IF160plus IF160mplus	IF260plus IF260mplus	IF450plus IF450mplus	IF750plus IF750mplus

Options	30	55	75	110	160	260	450	750
Voltage 115 V, 50/60 Hz					X2			
Extended overtemperature protection by additionally integrated Pt100 sensor for independent temperature monitoring for models with SingleDISPLAY					A6			
Chamber modification for the application of reinforced perforated stainless steel shelves or stainless steel grids (bearing rails mounted in the working chamber) – includes replacement of 2 standard grids by 2 reinforced grids					-			K1
Interior lighting for observing the load					R0			
Interior socket (can only be ordered with limited temperature range – max. +70 °C) ampacity 230 V, 2.2 A, can be switched off with the On/Off switch, cannot be switched individually, moisture tight IP68 (only with SingleDISPLAY), (option A8 necessary)					R3			
Entry port, 23 mm clear diameter, for introducing connections at the side, can be closed by flap and silicone stopper, standard positions	left centre/centre			F0				
	left centre/top			F1				
	right centre/centre			F2				
	right centre/top			F3				
Entry port, 23 mm clear diameter, can be closed by flap, in special positions (please state location)	left			F4				
	right			F5				
	rear			F6				
Entry port, 14 mm clear diameter, can be closed by flap, in special positions in the back wall (please, state location)					D6			
Entry port, 38 mm clear diameter, can be closed by flap, in special positions at the back (please, state location)					F7			
Entry port, 57 mm clear diameter, can be closed by flap, in special positions at the back (please, state location)					F8			
Entry port, 100 mm clear diameter, can be closed by flap, in special positions at the back (please, state location)	-		F9					
4 – 20 mA current loop interface (0 to +90 °C = 4 – 20 mA) Temperature controller, actual value Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring (max. 1 SingleDISPLAY, max. 3 TwinDISPLAY) – price per sensor					V3			
					V6			
Fan speed monitoring with switching-off the heating and with alarm in case of failure – only for IFplus/IFmplus					V4			
Works calibration certificate for 3 temperatures: +37 °C, +52 °C, +70 °C					D00126			
Door with lock (safety lock)					B6			
Door hinged on the left					B8			-
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28 for external monitoring (indicates when setpoint is reached)					H5			
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)					H6			
Potential-free contact (24 V/2 A) with socket to NAMUR NE 28, for signal generation, controlled by programme segment, for free-selectable functions to be activated (e.g. activation of audible and visual signals, exhaust motors, fans, stirrers, etc.). Only for units with TwinDISPLAY 2 contacts					H72			
Process-dependent programmable door lock (only for units with TwinDISPLAY)					D4			
Door-open-recognition, incl. alarm, shuts down fan and after 30 sec. also heating (only for units with TwinDISPLAY)					V5			
Flexible Pt100 for positioning in chamber or in load with socket, 4-pin, according to NAMUR NE 28, for external temperature recording (load temperature) max. 3 sensors					H4			
Flexible Pt100 temperature sensor, positioned flexibly in chamber or load, for local temperature measurement (up to 3 additional sensors are possible). The measured temperature can, if required, be indicated on the display, recorded in the integral data store, and can be documented via the AtmoCONTROL software					H8			
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6					C3			
Temperature restriction; Temperatures: +60, +70, +80, +95, +100, +120, +160, +180, +200 or +220 °C (Please, indicate upon ordering)					A8			
Castor frame (2-part), height 140 mm					R9			-

Accessories	30	55	75	110	160	260	450	750
Stainless steel grid, electropolished	E28884	E20164		E20165		E28891	E20182	
Additional reinforced stainless steel grid, electropolished, max. loading 60 kg; from size 450 with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber		-		E29767		E29766	B32190	
Perforated stainless steel shelf	B29727	B03916		B00325		B29725	B00328	
Additional reinforced stainless steel shelf, max. loading 60 kg; with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber		-					B32191	
Stainless steel slide-in drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	E02070	E02072		E02073		E29726	E02075	
Stainless steel slide-in drip tray, 15 mm rim, with guide bars and fixing screws (can be used only in connection with option K1)		-					B32763	
Stainless steel bottom drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	B04356	B04358		B04359		B29722	B04362	
Stainless steel bottom drip tray, 15 mm rim (can be used only in connection with option K1)		-					B34055	

Accessories	30	55	75	110	160	260	450	750
Wall bracket for wall mounting	B29755	B29756	B29757	B29758	B29759	-		
Guarantee extension by 1 year	GA1Q5					GA2Q5		
USB-Ethernet adapter	E06192							
Ethernet connection cable 5 m for computer interface	E06189							
USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation by unauthorised third parties. When reordering please specify serial number (only for units with TwinDISPLAY)	B33170							
USB stick with documentation software AtmoCONTROL and operation manual for products with SingleDISPLAY (the standard equipment of appliances with TwinDISPLAY includes one USB stick with AtmoCONTROL). When reordering please specify serial number	B33172							
Set of height adjustable feet (4 pcs)	B29768						-	
Stacking set (4 pcs) for stacking of appliances of same size	B29744					-		
Plug-in tube extension (outer diam. 60.3 mm, inner 57 mm), straight, for exhaust air ducting (if necessary for connection by hose)	B29718							
Plug-in tube extension (outer diam. 60.3 mm, inner 57 mm), angled, for exhaust air ducting (if necessary for connection by hose)	B29719							
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), with air slots	B29728	B29730	B29732	B29734	B29736	B29738	B29740	B29742
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), without air slots	B29729	B29731	B29733	B29735	B29737	B29739	B29741	B29743
Subframe, adjustable in height (size 30 to 75: height 600 mm, size 110 to 450: height 500 mm)	B29745	B29747		B29749		B29751	B29753	-
Subframe, on castors (size 30 to 75: height 660 mm, size 110 to 160: height 560 mm)	B29746	B29748		B29750		-		
Subframe, adjustable in height, height 130 mm, for example for units with fresh air filter	B33657	B33659		B33661		B33664	-	
Software conforming to FDA AtmoCONTROL. Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit (only for units with TwinDISPLAY). Respective IQ/OQ documents available in German and English language (without surcharge)	FDAQ1							
Integration of additional units (up to max. 15 units) into an already existent FDA-software licence (only for units with TwinDISPLAY)	FDAQ2							
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer	D00124							
IQ/OQ document with device-specific works test data for one free-selectable temperature value, incl. temperature distribution survey at Memmert for 9 measuring points (size 30), 27 measuring points (sizes 55 – 1060) to DIN 12880:2007-05. PQ check list as support for validation by customer. Price for further temperature values and validation at customer site on demand	D00125	D00127						



CO₂ Incubator ICOMed with TwinDISPLAY
Software AtmoCONTROL

Model sizes: 50 / 105 / 150 / 240

+18 °C to +50 °C

Humidity 40 to 97 % rh


CO₂ concentration 0 to 20 %

O₂ concentration 1 to 20 %

CO₂ INCUBATOR ICOMed Safety at all times. When it comes to safety and user friendliness, the highly modern CO₂ incubator ICOMed is the perfect solution: Thanks to the battery-buffered ControlCOCKPIT, the operating display, logging and CO₂ control remain fully functional even when there is a power failure. All parameters are logged in accordance with the FDA and, when individually adjusted ranges for CO₂, O₂, temperature and humidity are exceeded, notifications can be sent to a mobile phone in addition to an alarm.

The control technology is so finely tuned that the setpoint temperature is guaranteed to be reached without temperature overshoots. With its rounded corners, the interior is easy to clean and can be sterilised for 60 minutes at 180 °C (including all sensors).

All ICOMed models are classified as medical product class IIa.

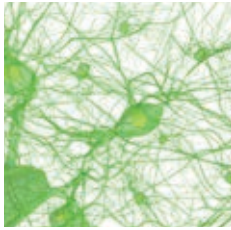

Unrivalled user friendliness

All parameters can be set easily and intuitively both with the ControlCOCKPIT or the AtmoCONTROL software. The shutter box can be opened, allowing fast access to controls. Maintenance is possible even if the appliances are stacked. The appliance has USB and Ethernet connections as well as a data logger with a ten-year storage capacity. Data can be read and programmes can be transferred by remote access.



Minimising vaporisation and condensation

The active humidity control minimises vaporisation in the interior and ensures short recovery times after the door has been opened. Together with the heating of the interior from all six sides including the heated inner glass door, it prevents the dangerous formation of condensation and offers maximum protection for cell and tissue cultures. The turbulence-free chamber ventilation ensures a constant and uniform atmosphere.



IVF module for models ICO50med/ICO105med

In order to keep vaporisation, condensation and recovery times at a minimum during in vitro fertilisation, the Petri dishes are cultivated in separate slide-in units. The slide-in units in the optional IVF module can be pulled out easily and with low vibration and are equipped with a pull-out lock.



The CO₂ Incubator ICOMed is a medical device:

Memmert subjected its CO₂ incubator ICOMed to a comprehensive medical device evaluation. Every Memmert CO₂ incubator ICOMed is classified as a Class IIa medical device. The ICOMed is intended for the creation and maintenance of constant environmental conditions for application in the field of in vitro fertilisation (IVF), especially for the incubation of oocytes, spermatozoa and zygotes in special culture dishes for IVF application as well as for gene expression and the biosynthesis of RNA and proteins. The CE label on the appliances includes the mark 0197, denoting TRLP – TÜV Rheinland as the notified body.




CO₂ INCUBATORS IC0med

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:



- Interior:** Stainless steel, material 1.4301 (ASTM 304), deep-drawn, seamlessly welded
- Housing:** Textured stainless steel, rear zinc-plated steel, intuitively operated TwinDISPLAY (TFT colour display) with touchscreen; fully insulated stainless steel door and heated inner glass door
- Automatic sterilisation:** Humidity and CO₂ sensor sterilised inside the CO₂ incubator
- Connection:** Mains cable with plug (German type)
- Installation:** 4 adjustable feet
- Interfaces:**

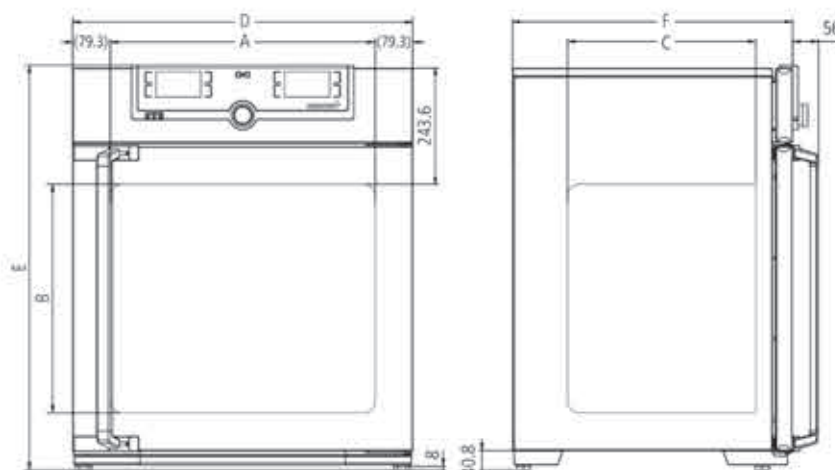

Model sizes/Description			50	105	150	240
Stainless steel interior	Volume	approx. l	56	107	156	241
	Width	(A) mm	400	560	600	600
	Height	(B) mm	425	480	700	810
	Depth (less 35 mm for fan)	(C) mm	330	400	500	500
	Max. number of perforated shelves	number	5	6	10	12
	Max. loading per perforated shelf	kg		15		
	Max. loading of chamber	kg	75	90	120	140
Textured stainless steel exterior	Width	(D) mm	559	719	759	759
	Height (variable through adjustable feet)	(E) mm	795	850	1070	1180
	Depth (without door handle, depth of door handle 56 mm)	(F) mm	521	591	691	691
	Fully insulated heated stainless steel door			●		
	Additional heated inner glass door			●		
Standard equipment	Stainless steel shelves, perforated	number	1	2		
	Stainless steel water dish (not applicable with option K7)			1		
	Works calibration certificate (measuring point chamber centre) at +37 °C, 5 % CO ₂ for standard units			●		
	Works calibration certificate at 37 °C, 5 % CO ₂ , 90 % rh and 10 % O ₂ (requires option K7 and option T6); standard equipment for units with O ₂ control			●		
	Works calibration certificate at 37 °C, 5 % CO ₂ and 90 % rh (requires option K7); standard equipment for units with active humidity control			●		
	CO ₂ connection set: hose with coupling and clamp			●		
	Standard sterilisation programme (without removing the sensors), humidity and CO ₂ sensor sterilised inside the CO ₂ incubator			60 minutes at 180 °C		
Temperature	Working temperature range	°C	5 above ambient temperature up to +50			
	Setting temperature range	°C	+18 to +50			
	Setting accuracy	°C	0.1			
	Temperature fluctuations with time at 37 °C (to DIN 12880:2007-05)	K	+/- 0.1			
	Temperature variation in chamber at +37 °C (to DIN 12880:2007-05)	K	+/- 0.3			
Humidity	Humidity limitation thanks to a Peltier element; when water dish is full and inserted, the Peltier element limits the value of relative humidity in the interior to 93 % rh +/- 2.5 %			●		
	Setting range active humidity control (with option K7)	% rh	40 to 97 and rh-Off			
	Setting accuracy	% rh	0.5			
CO ₂ / O ₂	Digital electronic CO ₂ control with dual beam NDIR system, with auto-diagnostic system and acoustic fault indication, barometric pressure compensation			●		
	Setting range CO ₂	% CO ₂	0 to 20			
	Variation in time CO ₂	% CO ₂	+/- 0.2			
	Setting accuracy CO ₂	% CO ₂	0.1			
	Setting range O ₂	% O ₂	1 to 20			
	Setting accuracy O ₂	% O ₂	0.1			
Further data	Electrical load at 230/115 V, 50/60 Hz	approx. W	1000	1500	2000	
Packing data	Net weight	approx. kg	55	75	90	110
	Gross weight (packed in carton)	approx. kg	74	100	116	145
	Width	approx. mm	730	800	840	840
	Height	approx. mm	950	1030	1250	1360
	Depth	approx. mm	640	800	900	900

Model sizes/Description	50	105	150	240
Order No. CO₂ Incubators	IC050med	IC0105med	IC0150med	IC0240med
Options	50	105	150	240
Voltage 115 V, 50/60 Hz		X2		
Battery-buffered ControlCOCKPIT: Uninterrupted supply for the entire display unit (ControlCOCKPIT) and therefore complete documentation of all parameters even when there is a power failure. The CO ₂ parameter is continuously regulated.		C2		
Two gas connections with quick release connectors for automatic switch-over of gas cylinders; incl. two CO ₂ connection sets: hose with coupling and clamp		T1		
Electropolished interior		T2		
Active microprocessor control for humidifying and dehumidifying (40 – 97 % rh), incl. digital indication and auto-diagnostic system ensures even more rapid reaching of set humidity and very short recovery times while avoiding condensate formation. Humidity supply with distilled water (from an external tank) by a self-priming pump; integral bacteria block by generating hotsteam, dehumidifying via sterile filter		K7		
Control of oxygen concentration by N ₂ inlet; adjustment range 1 % up to 20 % O ₂ ; setting accuracy 0.1 % (requires option K7). Incl. N ₂ connection set: hose with coupling and clamp		T6		
Peltier cooling unit enables a working temperature of 37 °C even at higher ambient temperatures of up to 35 °C		K5		
Capacitive humidity sensor for measuring and displaying the relative humidity		K6		
Entry port (silicone), 40 mm clear diameter, moisture tight, can be closed by silicone stopper, at the back, centre right; not available for IC050med with active humidity control (option K7) or humidity display (option K6)		F7		
Inner door with partitioned glass doors	-		K4	
4 – 20 mA current loop interface	Temperature controller, actual value (0 to +70 °C = 4 – 20 mA)	V3		
	Humidity controller, actual value (0 to 100 % rh = 4 – 20 mA) (requires option K7 or K6)	V7		
	CO ₂ controller, actual value (0 to 25 % CO ₂ = 4 – 20 mA)	V9		
	O ₂ controller, actual value (0 to 25 % O ₂ = 4 – 20 mA) (requires option T6)	V1		
Works calibration certificate for 5 %, 7 % and 10 % CO ₂ (measured at +37 °C) special works calibration certificates upon request		D00106		
Start-up of IC0Med incubators and brief training (D, A, CH only), through Memmert service, not subject to discount		K9		
Door hinged on the left		B8		
Potential-free contact (24 V/2 A) with socket to NAMUR NE 28 for external monitoring (indicates when set points of temperature and CO ₂ are reached)		H5		
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)		H6		
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6		C3		
MobileALERT for up to 4 alarm notifications; temperature and CO ₂ alarm (standard), additionally humidity alarm (when equipped with option K7) and O ₂ alarm (when equipped with option T6)		C4		
Accessories	50	105	150	240
Perforated stainless steel shelf	E35160	E37418		E35158
Water dish	B38737	B38000		
Subframe (622 mm high) adjustable in height (sizes 150/240: should not be used for 2 stacked units)	B33504	B33505		B33506
Subframe (130 mm high); sizes 150/240: only in combination with the corresponding stacking sets for stacked appliances	B33507	B33508		B33509
HEPA-filter for chamber (filter class E11) according to EN 1822, packed in sterile condition, incl. fixing unit		B38739		
CO ₂ pressure reducing valve to DIN 8546, incl. gas cylinder monitor		E02087		
N ₂ pressure reducing valve to DIN EN ISO 2503, incl. gas cylinder monitor (requires option T6)		E06162		
Central water supply, with filter cartridges for connection to the domestic water supply, only in combination with option K7. Product information on demand		ZWVR6		
Central water supply, without filter cartridges for connection to the domestic water supply (only for demineralised water in accordance with VDE 0510/DIN EN 50272), only in combination with option K7. Product information on demand		ZWVR7		
Guarantee extension by 1 year		GA3Q5		
Celltron benchtop shaker (not subject to discount) – accessories upon request	-	E06724		
IVF-module for IC050med: patented consisting of 6 slide-in units, a total of 12 special racks with indentations for 12 Petri dishes (60 mm diam.) resp. 24 Petri dishes (35 mm diam.), 2 racks with indentations for 3 special media tubes each; racks with indentations for 4-well dishes on demand; only for IC050med with the options K7 and F7; works calibration certificate (measuring point chamber centre) at +37 °C, 5 %, 6 % and 7 % CO ₂ as well as 90 % rh; 5 % O ₂ for IVF unit equipped with option T6	B44128	-		
IVF-module for IC0105med: patented, consisting of 8 slide-in units, a total of 16 special racks with indentations for 16 Petri dishes (60 mm diam.) resp. 32 Petri dishes (35 mm diam.), 2 racks with indentations for 3 special media tubes each; racks with indentations for 4-well dishes on demand; only for IC0105med with the options K7 and F7; works calibration certificate (measuring point chamber centre) at 37 °C, 5 %, 6 % and 7 % CO ₂ as well as 90 % rh; 5 % O ₂ for IVF unit equipped with option T6	-	B42398	-	
Holder for Petri dishes round (only in combination with IVF-module)	E37026		-	
Holder for Petri dishes square (only in combination with IVF-module)	E37308		-	
Holder for test tube (only in combination with IVF-module)	E37069		-	
Magnetic foil, can be labelled with a non-permanent board marker (only in combination with IVF-module)	E36651		-	
USB-Ethernet adapter		E06192		
Ethernet connection cable 5 m for computer interface		E06189		
USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation by unauthorised third parties. When reordering please specify serial number		B33170		
Stacking set (4 pcs) for stacking of appliances of same size	B29744		-	
Stacking set (consisting of stacking corners, one connecting plate for the rear, two wall brackets) for stacking of two units of same size	-	B42114	B42115	
Software conforming to FDA AtmoCONTROL. Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit. Respective IQ/OQ documents available in German and English language (without surcharge)		FDAQ1		
Integration of additional units (up to max. 15 units) into an already existent FDA-software licence		FDAQ2		

Accessories	50	105	150	240
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer				D00124
IQ/OQ document with device-specific works test data for one free-selectable CO ₂ , humidity and temperature value, incl. temperature distribution survey at Memmert for 27 measuring points to DIN 12880:2007-05, PQ check list as support for validation by customer (a free-selectable humidity value is only possible with option K7). Price for validation at customer site on demand				D38897
IQ/OQ document with device-specific works test data for one free-selectable CO ₂ and temperature value, incl. temperature distribution survey at Memmert for 27 measuring points to DIN 12880:2007-05, PQ check list as support for validation by customer. Price for validation at customer site on demand				D38898
External measuring instrument with additional measuring head for temperature and humidity measurement. Product information on demand				B04714



CO₂-cooled incubator ICPeco
with TwinDISPLAY
AtmoCONTROL software

Model sizes ICPeco: 260 / 450 / 750

-12 °C to +60 °C

Model size ICP: 55

0 °C to +60 °C

Model sizes ICP: 110 / 260 / 450 / 750

-12 °C to +60 °C

COMPRESSOR-COOLED INCUBATOR ICPeco These environmentally-friendly cooled incubators are cooled with climate-neutral CO₂. Thanks to this refrigerant's excellent thermodynamic properties and the finely adjusted control technology, an ICPeco is both powerful and high-precision. Without critical temperature overshoots, it keeps the temperatures exactly at the setpoint.




Refrigerant CO₂ is climate-neutral

A CO₂-cooled incubator ICP^{eco} is in every respect positive for the ecological balance of a laboratory. Legal restrictions for use are completely excluded in the future, as the refrigerant CO₂ (R744), unlike fluorine-based refrigerants, has no greenhouse gas reduction potential. It is a by-product of industrial processes, which is why far less energy is used for its manufacture than for synthetic, fluorinated refrigerants. R744 is neither flammable nor toxic, does not cause ozone depletion in the atmosphere and does not require disposal or recycling.

Refrigerant CO₂ ensures better cooling performance

The contribution to process optimisation is also impressive. An ICP^{eco} is virtually maintenance-free and extremely powerful. Compared to appliances with R134a as refrigerant, it has faster temperature change rates during cooling-down. Memmert cooled incubators ICP with refrigerant R134a will be available in parallel for a transitional period.

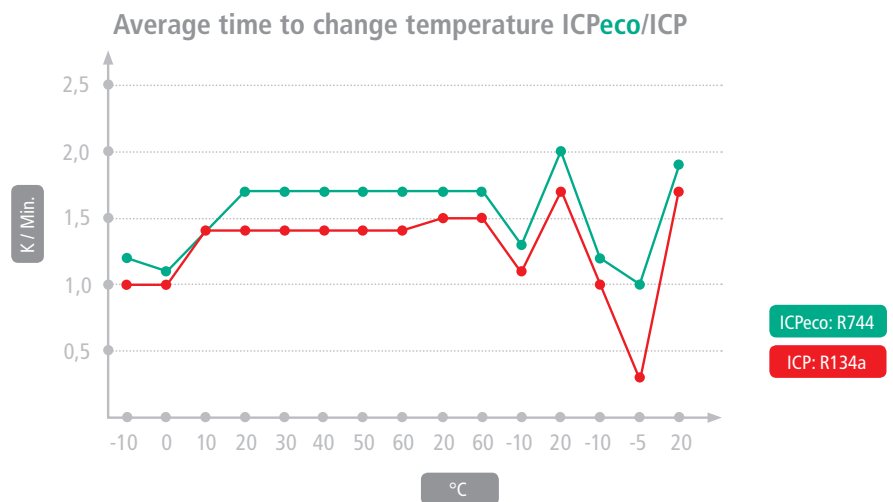


On average 20 % faster temperature change rates with a CO₂-cooled compressor (measurement ICP260^{eco} at ambient temperature +22 °C according to IEC 60068-3-5)





Completely enclosed working chamber

Cooling and heating units are situated outside the working chamber inside the air jacket temperature control system surrounding the entire chamber interior ensuring quick and precise temperature control. The motor-driven forced air circulation, adjustable in 10 % steps via the ControlCOCKPIT ensures optimum temperature distribution.

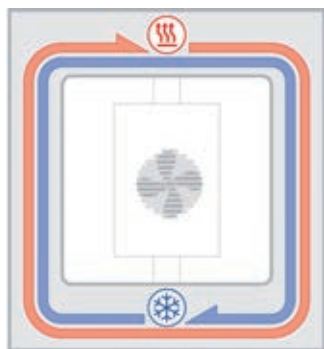

ICP air jacket temperature control system

Integrated energy saving function

The cooling unit works extremely energy-efficient because the heating is completely switched off in cooling mode. An intelligent DEFROST function enables defrosting as required.


COMPRESSOR-COOLED INCUBATORS ICPeco

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:



- Interior: Stainless steel, material 1.4301 (ASTM 304)
- Housing: Textured stainless steel, rear zinc-plated steel, intuitively operated TwinDISPLAY (TFT colour display) with touchscreen; inside glass door, outside fully insulated stainless steel door (from size 450 two leaves)
- Connection: Mains cable with plug (German type)
- Installation: Mounted on lockable castors
- Interfaces:

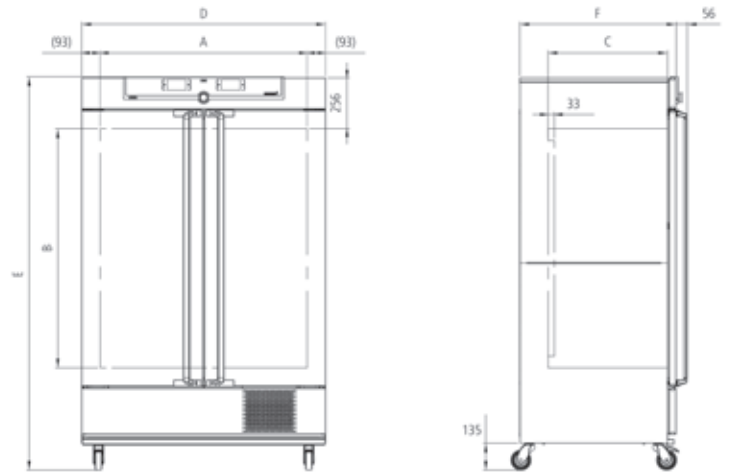

Model sizes/Description				260	450	750
Stainless steel interior	Volume		approx. l	256	449	749
	Width	(A)	mm	640	1040	
	Height	(B)	mm	800	720	1200
	Depth (less 33 mm for fan)	(C)	mm	500	600	
	Max. number of grids/shelves		number	9	8	14
	Max. loading per grid/shelf		kg	20	30	
	Max. loading of chamber		kg	200		
	Max. loading per slide-in drip tray		kg	4	8	
	Max. loading per bottom drip tray		kg	4	8	
Textured stainless steel exterior	Width	(D)	mm	824	1224	
	Height (with castors)	(E)	mm	1552	1613	1950
	Depth (without door handle), door handle + 56 mm	(F)	mm	684	784	
Standard equipment	Stainless steel grids, electropolished		number	2		
	Standard works calibration certificate (measuring point chamber center)		°C	+10 and +37		
Temperature	Working temperature range (not suitable for long-term storing at sub-zero temperatures. During permanent operation, the glass door may ice over)		°C	-12 to +60		
	Setting temperature range		°C	-12 to +60		
	Setting accuracy		°C	0.1		
Further data	Electrical load at 230 V, 50 Hz		approx. W	1200		
Packing data	Net weight		approx. kg	162	222	254
	Gross weight (packed in carton)		approx. kg	219	287	324
	Width		approx. mm	930	1330	
	Height		approx. mm	1760	1700	2150
	Depth		approx. mm	930	1050	

Order No. Compressor-Cooled Incubators

ICP260eco ICP450eco ICP750eco

Options	260	450	750
Chamber modification for the application of reinforced perforated stainless steel shelves or stainless steel grids (bearing rails mounted in the working chamber) – includes replacement of 2 standard grids by 2 reinforced grids	-	K1	
Interior socket, ampacity 230 V/2.2 A, can be switched off with the On/Off switch, cannot be switched individually, moisture tight IP68		R3	
Entry port, 23 mm clear diameter, for introducing connections at the side, can be closed by flap and silicone stopper, standard positions	left centre/centre	F0	
	left centre/top	F1	
	right centre/top	F3	
Entry port (silicone), 40 mm clear diameter, moisture tight, can be closed by silicone stopper, at the back (please, state location)		F7	
4 – 20 mA current loop interface	Temperature controller, actual value (-20 to +70 °C = 4 – 20 mA)	V3	
	Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring (max. 3) – price per sensor (-20 to +70 °C = 4 – 20 mA)	V6	
Fan speed monitoring with switching off the heating and with alarm in case of failure		V4	
Works calibration certificate for 3 temperatures: 0 °C, +37 °C, +60 °C		D00130	
Door with lock (safety lock)		B6	
Door hinged on the left	B8	-	
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28 for external monitoring (indicates when setpoint is reached)		H5	
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)		H6	
Potential-free contact (24 V/2 A) with socket to NAMUR NE 28, for signal generation, controlled by programme segment, for free-selectable functions to be activated (e.g. activation of audible and visual signals, exhaust motors, fans, stirrers, etc.)	2 contacts	H72	
Process-dependent programmable door lock		D4	
Door-open-recognition, incl. alarm		V5	
Flexible Pt100 for positioning in chamber or in load with socket, 4-pin, according to NAMUR NE 28, for external temperature recording (load temperature) max. 3 sensors		H4	
Flexible Pt100 temperature sensor, positioned flexibly in chamber or load, for local temperature measurement (up to 3 additional sensors are possible). The measured temperature can, if required, be indicated on the display, recorded in the integral data store, and can be documented via the AtmoCONTROL software		H8	
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6		C3	

Accessories	260	450	750
Stainless steel grid, electropolished	E28891	E20182	
Additional reinforced stainless steel grid, electropolished, max. loading 60 kg; from size 450 with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	E29766	B32190	
Perforated stainless steel shelf	B29725	B00328	
Additional reinforced stainless steel shelf, max. loading 60 kg; with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-	B32191	
Stainless steel slide-in drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	E29726	E02075	
Stainless steel slide-in drip tray, 15 mm rim, with guide bars and fixing screws (can be used only in connection with option K1)	-	B32763	
Stainless steel bottom drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	B29722	B04362	
Stainless steel bottom drip tray, 15 mm rim (can be used only in connection with option K1)	-	B34055	
USB-Ethernet adapter		E06192	
Ethernet connection cable 5 m for computer interface		E06189	
USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation by unauthorised third parties. When reordering please specify serial number		B33170	
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), with air slots	B29738	B29740	B29742
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), without air slots	B29739	B29741	B29743
Subframe, adjustable in height (height 500 mm)	B29751	B29753	-
Subframe, adjustable in height, height 130 mm, for example for units with fresh air filter	B33664	-	
Software conforming to FDA AtmoCONTROL. Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit. Respective IQ/OQ documents available in German and English language (without surcharge)		FDAQ1	
Integration of additional units (up to max. 15 units) into an already existent FDA-software licence		FDAQ2	
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer		D00124	
IQ/OQ document with device-specific works test data for one free-selectable temperature value, incl. temperature distribution survey at Memmert for 27 measuring points to DIN 12880:2007-05. PQ check list as support for validation by customer. Price for further temperature values and validation at customer site on demand		D00127	


COMPRESSOR-COOLED INCUBATORS ICP

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:



- Interior: Stainless steel, material 1.4301 (ASTM 304)
- Housing: Textured stainless steel, rear zinc-plated steel, intuitively operated TwinDISPLAY (TFT colour display) with touchscreen; inside glass door, outside fully insulated stainless steel door (from size 450 two leaves)
- Connection: Mains cable with plug (German type)
- Installation: Mounted on lockable castors
- Interfaces:

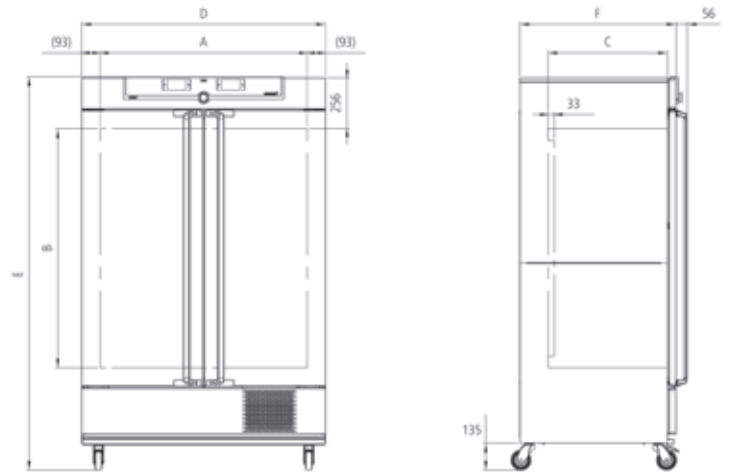

Model sizes/Description				55	110	260	450	750
Stainless steel interior	Volume	approx. l	53	108	256	449	749	
	Width	(A) mm	400	560	640	1040		
	Height	(B) mm	400	480	800	720	1200	
	Depth (less 33 mm for fan)	(C) mm	330	400	500	600		
	Max. number of grids/shelves	number	4	5	9	8	14	
	Max. loading per grid/shelf	kg	20			30		
	Max. loading of chamber	kg	80	150	200			
	Max. loading per slide-in drip tray	kg	1,5	3	4	8		
	Max. loading per bottom drip tray	kg	1,5	3	4	8		
Textured stainless steel exterior	Width	(D) mm	585	745	824	1224		
	Height (with castors)	(E) mm	1153	1233	1552	1613	1950	
	Depth (without door handle), door handle + 56 mm	(F) mm	514	584	684	784		
Standard equipment	Stainless steel grids, electropolished	number	1	2				
	Standard works calibration certificate (measuring point chamber center)	°C	+10 and +37					
Temperature	Working temperature range (not suitable for long-term storing at sub-zero temperatures. During permanent operation, the glass door may ice over)	°C	-12 to +60 (ICP55 0 to +60)					
	Setting temperature range	°C	-12 to +60 (ICP55 -5 to +60)					
	Setting accuracy	°C	0.1					
Further data	Electrical load at 230/115 V, 50/60 Hz	approx. W	1200					
Packing data	Net weight	approx. kg	89	113	157	217	249	
	Gross weight (packed in carton)	approx. kg	111	141	214	282	319	
	Width	approx. mm	760	880	930	1330		
	Height	approx. mm	1330	1410	1760	1700	2150	
	Depth	approx. mm	680	810	930	1050		

Order No. Compressor-Cooled Incubators

ICP55 ICP110 ICP260 ICP450 ICP750

Options	55	110	260	450	750
Voltage 115 V, 50/60 Hz	X2				
Chamber modification for the application of reinforced perforated stainless steel shelves or stainless steel grids (bearing rails mounted in the working chamber) – includes replacement of 2 standard grids by 2 reinforced grids	-			K1	
Interior socket, ampacity 230 V/2.2 A, can be switched off with the On/Off switch, cannot be switched individually, moisture tight IP68	R3				
Entry port, 23 mm clear diameter, for introducing connections at the side, can be closed by flap and silicone stopper, standard positions	left centre/centre		F0		
	left centre/top		F1		
	right centre/top		- F3		
Entry port (silicone), 40 mm clear diameter, moisture tight, can be closed by silicone stopper, at the back (please, state location)	F7				
4 – 20 mA current loop interface	4 – 20 mA current loop interface (-20 to +70 °C = 4 to 20 mA)		V3		
	Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring (max. 3) – price per sensor (-20 to +70 °C = 4 – 20 mA)		V6		
Fan speed monitoring with switching off the heating and with alarm in case of failure	V4				
Works calibration certificate for 3 temperatures: 0 °C, +37 °C, +60 °C	D00130				
Door with lock (safety lock)	B6				
Door hinged on the left	B8			-	
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28 for external monitoring (indicates when setpoint is reached)	H5				
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)	H6				
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28, triple, for signal generation, controlled by programme segment for a total of 3 freely selected functions to be activated (e.g. acoustic and visual signals, exhaust motors, fans, stirrers etc.).	2 contacts		H72		
Process-dependent programmable door lock	D4				
Door-open-recognition, incl. alarm	V5				
Flexible Pt100 for positioning in chamber or in load with socket, 4-pin, according to NAMUR NE 28, for external temperature recording (load temperature) max. 3 sensors	H4				
Flexible Pt100 temperature sensor, positioned flexibly in chamber or load, for local temperature measurement (up to 3 additional sensors are possible). The measured temperature can, if required, be indicated on the display, recorded in the integral data store, and can be documented via the AtmoCONTROL software	H8				
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6	C3				


Accessories	55	110	260	450	750
Stainless steel grid, electropolished	E20164	E20165	E28891	E20182	
Additional reinforced stainless steel grid, electropolished, max. loading 60 kg; from size 450 with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-	E29767	E29766	B32190	
Perforated stainless steel shelf	B03916	B00325	B29725	B00328	
Additional reinforced stainless steel shelf, max. loading 60 kg; with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-			B32191	
Stainless steel slide-in drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	E02072	E02073	E29726	E02075	
Stainless steel slide-in drip tray, 15 mm rim, with guide bars and fixing screws (can be used only in connection with option K1)	-			B32763	
Stainless steel bottom drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	B04358	B04359	B29722	B04362	
Stainless steel bottom drip tray, 15 mm rim (can be used only in connection with option K1)	-			B34055	
USB-Ethernet adapter			E06192		
Ethernet connection cable 5 m for computer interface			E06189		
USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation by unauthorised third parties. When reordering please specify serial number			B33170		
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), with air slots	B29730	B29734	B29738	B29740	B29742
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), without air slots	B29731	B29735	B29739	B29741	B29743
Subframe, adjustable in height (size 30 to 75: height 600 mm, size 110 to 450: height 500 mm)	B29747	B29749	B29751	B29753	-
Subframe, on castors (size 30 to 75: height 660 mm, size 110 to 160: height 560 mm)	B29748	B29750	-		
Subframe, adjustable in height, height 130 mm, for example for units with fresh air filter	B33659	B33661	B33664	-	
Software conforming to FDA AtmoCONTROL. Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit. Respective IQ/OQ documents available in German and English language (without surcharge)			FDAQ1		
Integration of additional units (up to max.15 units) into an already existent FDA-software licence (E05019)			FDAQ2		
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer			D00124		
IQ/OQ document with device-specific works test data for one free-selectable temperature value, incl. temperature distribution survey at Memmert for 27 measuring points to DIN 12880:2007-05. PQ check list as support for validation by customer. Price for further temperature values and validation at customer site on demand			D00127		



Peltier-cooled incubator IPP with SingleDISPLAY
Peltier-cooled incubator IPPplus with TwinDISPLAY
AtmoCONTROL software

Model sizes: 30 / 55 / 110 / 260 / 400 / 750 / 1060
0 °C to +70 °C

PELTIER-COOLED INCUBATOR IPP Heating and cooling seamlessly with one system thanks to Peltier technology. In this respect, cooled incubators IPP not only contribute to climate protection, but it also achieves an additional decrease in operating costs of up 90 % compared to compressor technology. This perfect development from the environmentally friendly and energy-saving heating/cooling technology by Memmert convinces by outstanding control precision and extremely small fluctuations.

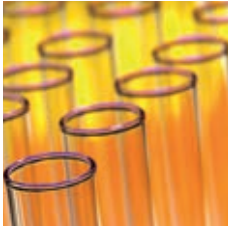

Extremely quiet and vibration-free

The fact that no compressor is required saves space and brings peace and quiet to the laboratory. As Peltier-cooled incubators IPP are almost vibration-free, they can also be applied in entomology. If defined humidity is also required, an alternative would be the constant climate chamber HPP, which is also equipped with Peltier technology.

No condensation in the interior chamber



Due to the closed Peltier cooling system, no outside air is exchanged. Physically derived, unavoidable formation of condensation during the cooling process does not take place in the interior chamber but on the outside heat sink. In addition, the in the Peltier elements integrated fans ensure a rapid transport of energy as well as an optimal temperature distribution.

Energy-saving heating/cooling technology combination

In contrast to compressor systems, Peltier technology is particularly economical at temperatures close to the ambient temperature, since energy is only required during heating or cooling. Therefore heating and cooling function are particularly precisely adjusted to each other.

Comparison compressor technology and Peltier technology

Reduction in energy consumption up to 90 %

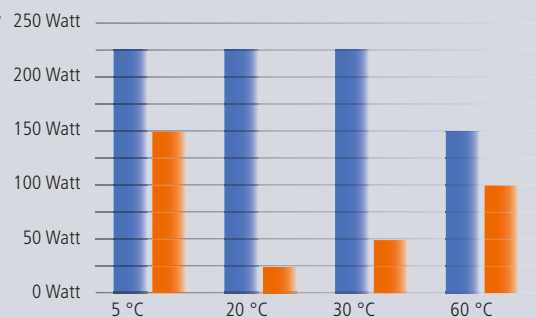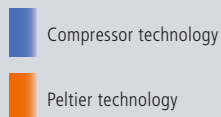

PELTIER-COOLED INCUBATORS IPP

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:
(EAC not valid for IPP400/1060)



Interior: Stainless steel, material 1.4301 (ASTM 304), deep-drawn

Housing: Textured stainless steel, rear zinc-plated steel, intuitively operated SingleDISPLAY or TwinDISPLAY (TFT colour display) with touchscreen

Double doors: Outside stainless steel, fully insulated, inside glass (size 750 and 1060 two leaves)


Connection: Mains cable with plug (German type)

Installation: 4 feet; sizes 400 to 1060 mounted on lockable castors

Interfaces:



USB: only for TwinDISPLAY

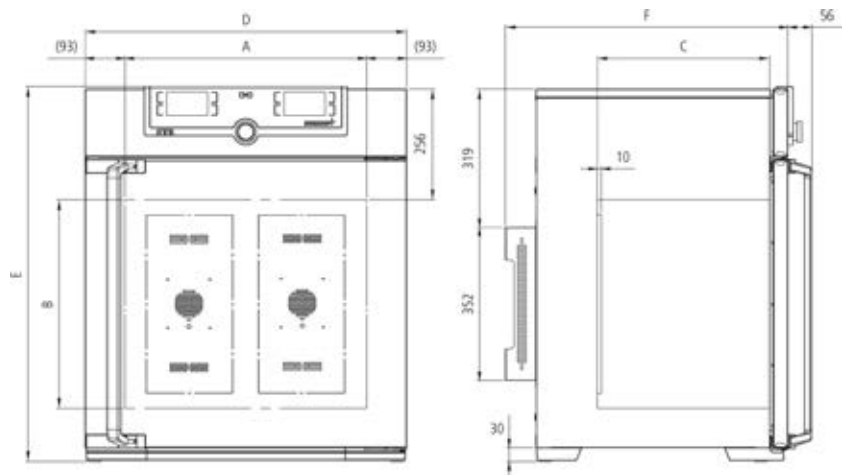

Model sizes/Description			30	55	110	260	400	750	1060
Stainless steel interior	Volume	approx. l	32	53	108	256	384	749	1060
	Width	(A) mm	400		560	640		1040	
	Height	(B) mm	320	400	480	800	1200		
	Depth (less 10 mm for fan – Peltier)	(C) mm	250	330	400	500		600	850
	Max. number of grids/shelves	number	3	4	5	9	14		
	Max. loading per grid/shelf	kg	20			30		20	
	Max. loading of chamber	kg	60	80	150	200			
	Max. loading per slide-in drip tray	kg	1,5		3	4	8		
	Max. loading per bottom drip tray	kg	1,5		3	4	8		
Textured stainless steel exterior	Width	(D) mm	585		745	824		1224	
	Height (sizes 400, 750, 1060 with castors)	(E) mm	704	784	864	1183	1720	1726	1661
	Depth (without door handle), door handle + 56 mm	(F) mm	506	586	656	756		856	1107
Standard equipment	Stainless steel grids, electropolished	number	1		2				
	Standard works calibration certificate (measuring point chamber center)	°C	+10 and +37						
Temperature	Working temperature range without light	°C	0 (at least 20 below ambient temperature) to +70						
	Working temperature range with light	°C	-		+10 to +40				-
	Setting temperature range	°C	0 to +70						
	Setting accuracy	°C	0.1						
Further data	Electrical load at 230/115 V, 50/60 Hz	approx. W	140	275	550	820	1100	1300	1500
	Peltier elements in the rear	number	1		2	3	5	6	
Packing data	Net weight	approx. kg	40	52	78	114	157	230	255
	Gross weight (packed in carton)	approx. kg	56	71	103	165	210	301	419
	Width	approx. mm	660	730	830	930		1330	1370
	Height	approx. mm	890	950	1050	1380	1930	1910	1970
	Depth	approx. mm	650	670	800	930		1050	1300

Order No. Peltier-Cooled Incubators

IPP = Peltier-Cooled Incubators

plus = Model with TwinDISPLAY

IPP30	IPP55	IPP110	IPP260	IPP400	IPP750	IPP1060
IPP30plus	IPP55plus	IPP110plus	IPP260plus	IPP400plus	IPP750plus	IPP1060plus

Options	30	55	110	260	400	750	1060
Voltage 115 V, 50/60 Hz	X2						
Chamber modification for the application of reinforced perforated stainless steel shelves or stainless steel grids (bearing rails mounted in the working chamber) – includes replacement of 2 standard grids by 2 reinforced grids	-					K1	-
Light module cold white 6,500 K: light strips arranged on the side walls of the interior, 10 strips for model 110, 14 for model 260/400/750, programme-controlled dimming from 0 to 100 % (in 1 % steps), ramp programming in combination with temperature (only with TwinDISPLAY)	-	T7					-
Light module cold white 6,500 K + warm white 2,700 K: LED light strips – 10 strips for model 110, 14 for model 260/400/750 – (5 resp. 7 alternating cold white light strips and 5 resp. 7 warm white light strips) on the side walls of the interior, programme-controlled dimming from 0 to 100 % (in 1 % steps), ramp programming in combination with temperature (only with TwinDISPLAY)	-	T8					-
Light module warm white 2,700 K: light strips arranged on the side walls of the interior, 10 strips for model 110, 14 for model 260/400/750, programme-controlled dimming from 0 to 100 % (in 1 % steps), ramp programming in combination with temperature (only with TwinDISPLAY)	-	T9					-
Interior socket, ampacity 230 V/2.2 A, can be switched off with the On/Off switch, cannot be switched individually, moisture tight IP68	R3						
Entry port, 23 mm clear diameter, for introducing connections at the side, can be closed by flap, standard positions (F0 and F2 not for model size 260 with light module; F0 – F3 not for model size 110 with light module)	left centre/centre left centre/top right centre/centre right centre/top		F0 F1 F2 F3				
Entry port, 23 mm clear diameter, can be closed by flap (please, state location)	left right rear		F4 F5 F6				
Entry port, 14 mm clear diameter, can be closed by flap, in special positions in the back wall (please, state location)	D6						
Entry port, 38 mm clear diameter, can be closed by flap, in special positions in the back wall (please, state location)	F7						
4 – 20 mA current loop interface (-10 to +80 °C = 4 – 20 mA)	Temperature controller, actual value Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring (max. 1 SingleDISPLAY, max. 3 TwinDISPLAY) – price per sensor			V3 V6			
Works calibration certificate for 3 temperatures: +5 °C, +37 °C, +60 °C	D00129						
Door with lock (safety lock)	B6						
Door hinged on the left	B8					-	
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28 for external monitoring (indicates when setpoint is reached)	H5						
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)	H6						
Potential-free contact (24 V/2 A) with socket to NAMUR NE 28, for signal generation, controlled by programme segment, for free-selectable functions to be activated (e.g. activation of audible and visual signals, exhaust motors, fans, stirrers, etc.). Only for units with TwinDISPLAY	2 contacts		H72				
Process-dependent programmable door lock (only for units with TwinDISPLAY)	D4						
Door-open-recognition, incl. alarm, shuts down fan and after 30 sec. also heating (only for units with TwinDISPLAY)	V5						
Flexible Pt100 for positioning in chamber or in load with socket, 4-pin, according to NAMUR NE 28, for external temperature recording (load temperature) max. 3 sensors	H4						
Flexible Pt100 temperature sensor, positioned flexibly in chamber or load, for local temperature measurement (up to 3 additional sensors are possible). The measured temperature can, if required, be indicated on the display, recorded in the integral data store, and can be documented via the AtmoCONTROL software	H8						
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6	C3						
Castor frame (2-part), height 140 mm	R9					-	

Accessories	30	55	110	260	400	750	1060
Stainless steel grid, electropolished	E28884	E20164	E20165	E28891	E20182	B41251	
Additional reinforced stainless steel grid, electropolished, max. loading 60 kg; size 750 with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-		E29767	E29766	B32190	-	
Perforated stainless steel shelf	B29727	B03916	B00325	B29725	B00328	B32549	
Additional reinforced stainless steel shelf, max. loading 60 kg; with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-					B32191	-
Stainless steel slide-in drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	E02070	E02072	E02073	E29726	E02075	B32599	
Stainless steel slide-in drip tray, 15 mm rim, with guide bars and fixing screws (can be used only in connection with option K1)	-					B32763	-
Stainless steel bottom drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	B04356	B04358	B04359	B29722	B04362	B29769	
Stainless steel bottom drip tray, 15 mm rim (can be used only in connection with option K1)	-					B34055	-
Guarantee extension by 1 year	GA1Q5			GA2Q5		GA4Q5	
USB-Ethernet adapter				E06192			
Ethernet connection cable 5 m for computer interface				E06189			
USB User-ID stick (with User-ID licence): Oven-linked authorisation licence (User-ID-programme) on Memory-stick, prevents undesired manipulation by unauthorised third parties. When reordering please specify serial number (only for units with TwinDISPLAY)				B33170			

Accessories	30	55	110	260	400	750	1060
USB stick with documentation software AtmoCONTROL and operation manual for products with SingleDISPLAY (the standard equipment of appliances with TwinDISPLAY includes one USB stick with AtmoCONTROL). When reordering please specify serial number	B33172						
Set of height adjustable feet (4 pcs)	B29768				-		
Stacking set (4 pcs) for stacking of appliances of same size	B29744			-			
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), with air slots	B29728	B29730	B29734	B29738	B42116	B29742	
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), without air slots	B29729	B29731	B29735	B29739	B42117	B29743	
Subframe, adjustable in height (size 30 to 75: height 600 mm, size 110 to 450: height 500 mm)	B29745	B29747	B29749	B29751	-		
Subframe, on castors (size 30 to 75: height 660 mm, size 110 to 160: height 560 mm)	B29746	B29748	B29750	-			
Subframe, adjustable in height, height 130 mm, for example for units with fresh air filter	B33657	B33659	B33661	B33664	-		
Software conforming to FDA AtmoCONTROL. Meets the requirements for the use of electronically stored data sets and electronic signatures as laid down in Regulation 21 CFR Part 11 of the US Food and Drug Administration (FDA). Base licence for the control of one unit (only for units with TwinDISPLAY). Respective IQ/OQ documents available in German and English language (without surcharge)	FDAQ1						
Integration of additional units (up to max. 15 units) into an already existent FDA-software licence (only for units with TwinDISPLAY)	FDAQ2						
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer	D00124						
IQ/OQ document with device-specific works test data for one free-selectable temperature value, incl. temperature distribution survey at Memmert for 9 measuring points (size 30), 27 measuring points (sizes 55 – 1060) to DIN 12880:2007-05. PQ check list as support for validation by customer. Price for further temperature values and validation at customer site on demand	D00125	D00127					
External measuring instrument with sensors for daylight and UV-light. Product information on demand (models IPPplus)	B04713						-
External measuring instrument with additional measuring head for temperature and humidity measurement. Product information on demand (models IPPplus)	B04714						-



Cooled storage incubator IPS
with SingleDISPLAY
AtmoCONTROL software

Model sizes: 260 / 750
+14 °C to +45 °C

COOLED STORAGE INCUBATOR IPS Save energy and reduce the strain on the climate at the same time! If microbiological cultures, BOB5 samples, drinks containers or cosmetics need to be stored over a long period at constant temperatures, cooled storage incubators IPS with energy-efficient Peltier technology are the perfect choice: absolute reliability, precision, durability and eco-friendliness.



Considerable potential for savings in acquisition and operating costs

Temperature changes are not always necessary for long-term storage or incubating. So why design heating, cooling and controlling systems for rapid heating up and cooling down times? The performance of the IPS was tailor-made for permanent operation at constant temperatures close to room temperature. The advantage: Acquisition costs and operating costs are considerably reduced in comparison to conventional cooled incubators with compressor technology, as well as to a large Peltier-cooled incubator.



Ideal for high ambient temperatures

Thanks to Peltier elements integrated for cooling the working chamber, the chamber load won't break into sweat even at high ambient temperatures. Constant and precise incubation at room temperature is guaranteed.



Low in vibration and durable for absolutely safe long-term storage

Like the cooled incubator IPP, the IPS offers all the advantages of Peltier technology to the user. Its interior chamber that is completely insulated from the environment minimises the risk of drying out of the samples. It is practically noise-free and not only reduces stress on the chamber load but also soothes the nerves of employees thanks to its quiet operation.



Glimpse into a Memmert storage incubator:
Peltier elements guarantee perfect climate inside the chamber.




COOLED STORAGE INCUBATORS IPS

according to DIN 12880:2007-05, EN 61010-1 (IEC 61010-1), EN 61010-2-010

Standard units are safety-approved and bear the test marks:



- Interior: Stainless steel, material 1.4301 (ASTM 304), deep-drawn
- Housing: Textured stainless steel, rear zinc-plated steel, intuitively operated SingleDISPLAY (TFT colour display) with touchscreen
- Double doors: Outside stainless steel, fully insulated, inside glass (size 750 two leaves)
- Connection: Mains cable with plug (German type)
- Installation: 4 feet; size 750 mounted on lockable castors
- Interfaces:

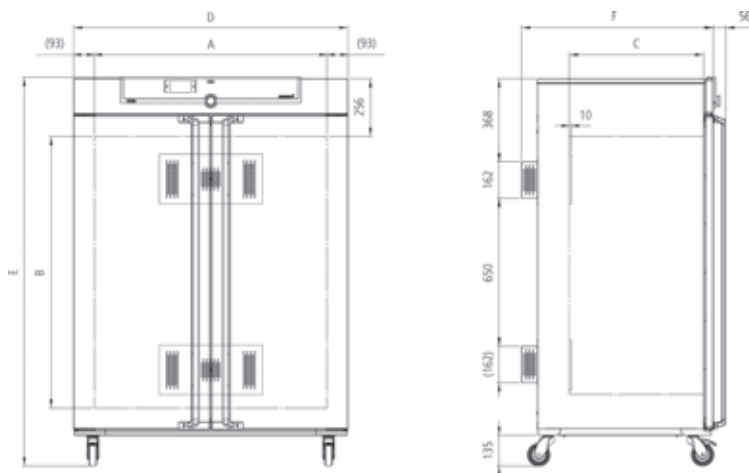

Model sizes/Description			260	750
Stainless steel interior	Volume	approx. l	256	749
	Width	(A) mm	640	1040
	Height	(B) mm	800	1200
	Depth (less 10 mm for fan – Peltier)	(C) mm	500	600
	Max. number of grids/shelves	number	9	14
	Max. loading per grid/shelf	kg	20	30
	Max. loading of chamber	kg	200	
	Max. loading per slide-in drip tray	kg	4	8
	Max. loading per bottom drip tray	kg	4	8
Textured stainless steel exterior	Width	(D) mm	824	1224
	Height (size 750 with castors)	(E) mm	1183	1726
	Depth (without door handle), door handle + 56 mm	(F) mm	754	856
Standard equipment	Stainless steel grids, electropolished	number	2	
	Standard works calibration certificate (measuring point chamber center)	°C	+18 and +25	
Temperature	Working temperature range	°C	+14 to +45	
	Setting temperature range	°C	+14 to +45	
	Setting accuracy	°C	0.1	
Further data	Electrical load at 230/115 V, 50/60 Hz	approx. W	550	
	Peltier elements in the rear	number	2	
Packing data	Net weight	approx. kg	113	230
	Gross weight (packed in carton)	approx. kg	164	301
	Width	approx. mm	930	1330
	Height	approx. mm	1380	1910
	Depth	approx. mm	930	1050
Order No. Cooled Storage Incubators			IPS260	IPS750

Options	260	750
Voltage 115 V, 50/60 Hz		X2
Chamber modification for the application of reinforced perforated stainless steel shelves or stainless steel grids (bearing rails mounted in the working chamber) – includes replacement of 2 standard grids by 2 reinforced grids	-	K1
Interior socket, ampacity 230 V/2.2 A, can be switched off with the On/Off switch, cannot be switched individually, moisture tight IP68		R3
Entry port, 23 mm clear diameter, for introducing connections at the side, can be closed by flap, standard positions	left centre/centre	F0
	left centre/top	F1
	right centre/centre	F2
	right centre/top	F3
Entry port, 23 mm clear diameter, can be closed by flap, in special positions (please state location)	left	F4
	right	F5
	rear	F6
Entry port, 14 mm clear diameter, can be closed by flap, in special positions in the back wall (please, state location)		D6
Entry port, 38 mm clear diameter, can be closed by flap, in special positions in the back wall (please, state location)		F7
4 – 20 mA current loop interface (0 to +70 °C = 4 – 20 mA)	Temperature controller, actual value	V3
	Temperature of a Pt100 sensor positioned flexibly in chamber for external temperature monitoring	V6
Works calibration certificate for a freely selectable temperature value		D00131
Door with lock (safety lock)		B6
Door hinged on the left	B8	-
Potential-free contact (24 V/2 A) with socket, according to NAMUR NE 28 for external monitoring (indicates when setpoint is reached)		H5
Potential-free contact for combination error message (e.g. supply failure, sensor fault, fuse)		H6
Flexible Pt100 for positioning in chamber or in load with socket, 4-pin, according to NAMUR NE 28, for external temperature recording (load temperature)		H4
Flexible Pt100 temperature sensor, positioned flexibly in chamber or load, for local temperature measurement (up to 3 additional sensors are possible). The measured temperature can, if required, be indicated on the display, recorded in the integral data store, and can be documented via the AtmoCONTROL software		H8
MobileALERT, notification by SMS in case of any error or alarm of the device. Requires option H6		C3
Castor frame (2-part), height 140 mm	R9	-

Accessories	260	750
Stainless steel grid, electropolished	E28891	E20182
Additional reinforced stainless steel grid, electropolished, max. loading 60 kg; size 750 with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	E29766	B32190
Perforated stainless steel shelf	B29725	B00328
Additional reinforced stainless steel shelf, max. loading 60 kg; with guide bars and fixing screws (only in connection with option K1). Please consider max. loading of chamber	-	B32191
Stainless steel slide-in drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	E29726	E02075
Stainless steel slide-in drip tray, 15 mm rim, with guide bars and fixing screws (can be used only in connection with option K1)	-	B32763
Stainless steel bottom drip tray, 15 mm rim (may affect the temperature distribution) – cannot be used in connection with option K1	B29722	B04362
Stainless steel bottom drip tray, 15 mm rim (can be used only in connection with option K1)	-	B34055
Guarantee extension by 1 year		GA2Q5
USB-Ethernet adapter		E06192
Ethernet connection cable 5 m for computer interface		E06189
USB stick with documentation software AtmoCONTROL and operation manual. When reordering please specify serial number.		B33172
Set of height adjustable feet (4 pcs)	B29768	-
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), with air slots	B29738	B29742
Flush-fit unit (stainless steel frame covering gap between oven and wall opening), without air slots	B29739	B29743
Subframe, adjustable in height (height 500 mm)	B29751	-
Subframe, adjustable in height, height 130 mm, for example for units with fresh air filter	B33664	-
IQ document with device-specific works test data, OQ/PQ check list as support for validation by customer		D00124
IQ/OQ document with device-specific works test data for one free-selectable temperature value, incl. temperature distribution survey at Memmert for 27 measuring points to DIN 12880:2007-05. PQ check list as support for validation by customer. Price for further temperature values and validation at customer site on demand		D00127

SOFTWARE AtmoCONTROL


AtmoCONTROL

The innovative control and logging software

Parameters such as temperature and humidity as well as the process time can be set directly at the ControlCOCKPIT. Ramp programming is done via the control and logging software AtmoCONTROL, which features a completely new software design.

Drag, drop & go!

Numerical and graphic programming of complex processes is a thing of the past. Today, programming is done via AtmoCONTROL by means of the mouse or touchpad on your notebook. Even the most complex ramp programmes are created within minutes. Simply drag & drop the graphical symbols for the desired parameters to the input field and change the values according to your wishes with a mouse click.

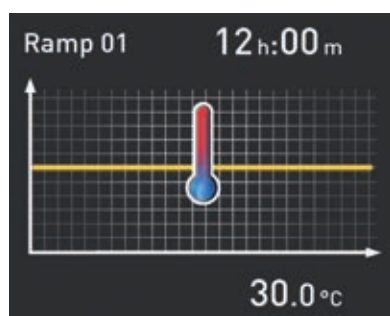

Programme functions for appliances with SingleDISPLAY and TwinDISPLAY

- Reading out, managing and organising the data logger
- Saving the log memory in various formats
- Online monitoring of up to 32 connected appliances
- Optical alarms when the alarm limits individually set at the ControlCOCKPIT are exceeded
- Automatic alarm to one or several e-mail addresses

Additional functions for appliances with TwinDISPLAY

- Intuitive programming and archiving of ramps and programme sequences
- Synchronous visualisation of the created programme sequence during programming
- Application-specific repeat functions (loops) can be inserted within a temperature control programme in any place
- Simple creation of repeating weekly programmes
- Programming, managing and transferring programmes via Ethernet interface or USB port



MODEL VARIANTS

SingleDISPLAY ControlCOCKPIT with one TFT display	TwinDISPLAY ControlCOCKPIT with two TFT displays
AVAILABLE APPLIANCES UN/UNm / UF/UFm / IN/INm / IF/IFm / IFbw / SN / SF / IPP / IPS	AVAILABLE APPLIANCES UNplus/UNmplus / UFplus/UFmplus / UF TS / UNpa INplus/INmplus / IFplus/IFmplus / SNplus / SFplus / VO ICOMed / IPPplus / ICPeco / ICP / HPP / ICHeco / ICH / HCP
One high-resolution TFT colour display with touch-sensitive buttons for selection of functions	Two high-resolution TFT colour displays with touch-sensitive buttons for selection of functions
Available parameters on the ControlCOCKPIT: Temperature (Celsius or Fahrenheit), fan speed, exhaust air flap position, programme time	Available parameters on the ControlCOCKPIT: Temperature (Celsius or Fahrenheit), fan speed, exhaust air flap position, programme time, relative humidity, illumination, CO ₂
One temperature sensor Pt100 DIN class A in a 4-wire circuit	Two Pt100 sensors DIN class A in a 4-wire circuit for mutual monitoring, taking over functions in case of an error
	HeatBALANCE function for application specific adjustment of heat output distribution (balance) between the upper and lower heating groups in an adjustment range between -50 % and +50 % (not valid for models 30, HPP110, IPP110plus, ICP, ICH)
AtmoCONTROL software for reading out, managing and organising the data logger via Ethernet interface (temporary trial version can be downloaded). USB stick with AtmoCONTROL software available as accessory (on demand)	AtmoCONTROL software on a USB stick for programming, managing and transferring programmes via Ethernet interface or USB port
	ControlCOCKPIT with USB port for uploading programmes, reading out protocol logs, activating the User-ID function
	Displaying of already logged protocol data on the ControlCOCKPIT (max 10,000 values correspond to approx. 1 week)
Ethernet interface on the rear of the appliance for reading out the protocol log and for online logging	Ethernet interface on the rear of the appliance for reading out the protocol log and for uploading programmes and for online logging
Double overtemperature protection: Electronic temperature monitoring with freely adjustable monitoring temperature, for models U, I, S with option A6 TWW/TWB (protection class 3.1 or 2), mechanical temperature limiter TB acc. to DIN 12880	Multiple overtemperature protection: Electronic temperature monitoring TWW/TWB (protection class 3.1 or 2 resp. 3.3 for units with active cooling) and mechanical temperature limiter TB (protection class 1) acc. to DIN 12880, AutoSAFETY automatically adjusts to the set value within a freely adjustable tolerance range. Setting individual MIN / MAX values for over/undertemperature alarm and also for all other parameters such as relative humidity, CO ₂
PID microprocessor control with integrated auto-diagnostic system	
Structured stainless steel housing, scratch-resistant, robust and durable; rear of zinc-plated steel	
High-temperature connectors on the rear of the appliance for single-phase power connection according to country specific systems and IEC standards	
Internal data logger with a storage capacity of at least 10 years	
German, English, French, Spanish, Polish, Czech, Hungarian language settings available on the ControlCOCKPIT	
Digital backwards counter with target time setting, adjustable from 1 minute to 99 days	
The SetpointWAIT function guarantees that the process time does not start until the set temperature is reached at all measuring points – optional for temperature values recorded by the freely positionable Pt100 sensors inside the chamber	
Adjustment of three calibration values for temperature and additional appliance specific parameters directly at the ControlCOCKPIT	



memmert
Experts in Thermostatics

HEATING AND DRYING OVENS

UNIVERSAL OVEN U

PASS-THROUGH OVEN UF TS

PARAFFIN OVEN UNpa

STERILISER S

VACUUM OVEN VO

BLANKET WARMER IFbw

INCUBATORS

INCUBATOR I

CO₂ INCUBATOR ICOfed

COMPRESSOR-COOLED INCUBATOR ICPecc/ICP

PELTIER-COOLED INCUBATOR IPP

COOLED STORAGE INCUBATOR IPS

CLIMATE CHAMBERS

CONSTANT CLIMATE CHAMBER HPP

HUMIDITY CHAMBER HCP

CLIMATE CHAMBER ICHeco/ICH

ENVIRONMENTAL TEST CHAMBER CTC/TTC

WATERBATHS / OILBATHS

WATERBATH W

OILBATH O

Distributed and Supported in the UK by:



24 Norman Way Industrial Estate, Over, Cambridge, CB24 5WE
Tel: +44(0)1954 233 100 Fax: +44(0)1954 233 101
Email: sales@camlab.co.uk Web: www.camlab.co.uk



Memmert GmbH + Co. KG
P.O. Box 1720 | D-91107 Schwabach
Tel. +49 9122 925-0 | Fax +49 9122 14585
E-Mail: sales@memmert.com
facebook.com/memmert.family
The platform for experts: www.atmosafe.net