

Product Information. **PRODUCT CODE: 1181400**

Big Blue Wipes are effective against 99.9% of gram-positive and gram-negative bacteria, as well as most viruses, fungi and other micro-organisms including spores. They are pre-dosed with a dual active formulation, providing effective infection control and preventing cross contamination of surfaces.

Areas of use:

This large bucket of 1000 disinfectant wipes has been designed for low cost disinfection in high use areas, ideal for the disinfection of large surfaces, machinery and equipment. Pal Disinfectant wipes can be used in potentially sensitive areas including food production, food service, healthcare, general industry and public environments.

Benefits of use:

Compliant

- Controlled dosage ensures effective disinfection for every use.
- Chemical solutions proven effective against a wide range of organisms.
- Can be used as part of a clean-down programme.

Convenient

- No requirement for pre-mixing, reducing wastage and risk of errors.
- Large bucket with carry handle for easy transportation.
- Wall brackets available for extra convenience.

Cost effective

- Saves valuable time and reduces the need for buckets and cloths.
- Refills are available bringing additional cost savings.
- Large bucket of 1000 wipes means low cost per wipe.

Use biocides safely. Always read the label and product information sheet before use.



Product Specification.

Wipe Material:

26gsm Thermo-bonded, Polypropylene and Viscose nonwoven.

Quantity/Tub: 1000 wipes

Outer case Quantity: N/A

Wipe Size: 195mm x 200mm

Material Colours:

Chemical Solution:

Surface Disinfectant

Ingredients:

Alkyl Dimethyl Benzyl, Ammonium Chloride
Alkyl Dimethyl Ethylbenzyl
Ammonium Chloride (Quaternary,
Ammonium Chlorides) 0.48%; Polymeric
Biguanide Hydrochloride 0.2%

Testing and Certifications:

Independently certified to kill a wide range of Bacteria, Viruses, Fungi and Spores – see attached technical sheet for details.

Shipping & Pallet Information.

Product weight: 3.4kg

Product dimensions: 270 x 270 x 230mm

Product cube: 0.016m

Layers: 6

Cases/layer: 13


Pallet quantity: 78

Pallet type: Euro

Pallet weight: 280kg

Dangerous Goods

Information: N/A

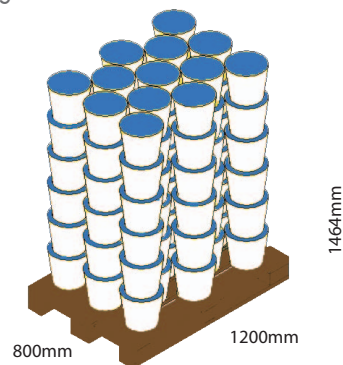

Pal Disinfectant wipes are developed and tested in our own laboratories by qualified chemists and microbiologists. They are then subjected to stringent EN testing procedures by independent testing laboratories to ensure efficacy against a wide range of Bacteria, Viruses, Fungi and Spores. Certificates are available upon request.

This table is a summary, listing the organisms tested and the EN Standards achieved. Where some organisms have been tested against more than one standard, results may have been omitted to avoid repetition.

			TEST CONDITIONS	
EN STANDARD TESTS	MICRO ORGANISMS	CONTACT TIME	CLEAN	DIRTY
BACTERIOCIDAL				
EN 1276 (Suspension Test) Log 5 kill in 5 minutes	Clostridium.difficile (vegetative)	5 mins	✓	
	Escherichia.coli	5 mins		✓
	Clostridium.perfringens (vegetative)	5 mins	✓	
	Listeria.monocytogenes	5 mins	✓	
	Proteus.vulgaris	1 min	✓	
	Salmonella.choleraesuis	5 mins	✓	
EN 13727 (Suspension Test) Log 5 kill in 60 minutes	Staphylococcus.aureus	1 min		✓
	Pseudomonas.aeruginosa	1 min		✓
	Enterococcus.hirae	1 min		✓
	MRSA	1 min		✓
	VRE	1 min		✓
EN 13697 (Surface Test) Log 4 kill in 5 minutes	Escherichia.coli	5 mins	✓	
	Campylobacter.jejuni	5 mins		✓
	Klebsiella.pneumoniae/NDM1	1 min		✓
EN 14561 (Surface Test) Log 5 kill in 60 minutes	Staphylococcus.aureus	5 mins		✓
	Pseudomonas.aeruginosa	5 mins		✓
	Enterococcus.hirae	5 mins		✓
TUBERCULOCIDAL				
EN 14348 (Suspension Test) Log 4 kill in 60 minutes	Mycobacterium.terrae	5 mins		✓
SPORICIDAL				
EN 13704 (Suspension Test) Log 3 kill in 60 minutes	Bacillus.subtilis spores	30 secs		✓
	Bacillus.cereus spores	30 secs		✓
VIRUCIDAL				
Blu Test Protocol Log 3 kill	Influenza A H1N1	5 mins	✓	
	BVDV (Hepatitis C surrogate)	5 mins	✓	
	FIV Petaluma (HIV surrogate)	5 mins	✓	
	Feline Coronavirus (SARS surrogate)	30 mins	✓	
YEASTICIDAL/FUNGICIDAL				
EN 14562 (Surface Test) Log 4 kill in 60 minutes	Candida albicans	15 mins		✓
	Aspergillus niger	15 mins		✓

Product has exceeded the required standard.