

## DIRECTIONS FOR USE: PORTABLE SHOWER STATION

### UTILIZATION:



1. Remove the safety pin
2. Press the operating lever
3. Press spraying trigger

### MAINTENANCE:




#### FILLING UP :

Unscrew 3 brass screw with a pin-spanner n°13 and remove the under-steam part.

Fill the water space with mineral water only  
Place the under-steam part and screw.

#### CHANGING CO<sub>2</sub> CARTRIDGE:

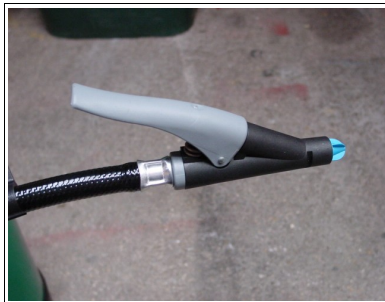

Check if the safety pin locks the operating lever.

Unscrew 3 brass screw with a pin-spanner n°13 and remove the under-steam part.

Unscrew the old CO<sub>2</sub> cartridge totally to remove it. Then place the new cartridge and screw it totally.

Place the under-steam part and screw

#### ATTENTION :

Before unscrewing, check if the shower is not under steam.

If it is under steam, press the safety pin to evacuate the whole CO<sub>2</sub>

**VERIFICATION:**



Gaz cartridge must be checked every year to ensure:

- There is no oxyde trace, which is dangerous, and can create a leakage.
- The CO<sub>2</sub> capacity is in the safety limit.

We recommend to contact a skilled expert to do this control once a year.

**ATTENTION:**

- Changing mineral water every 6 months.
- CO<sub>2</sub> cartridge must be changed after each utilization.