

CUB EASY GLASS WASHER PROGRAMS



Description of the phases in each program

Prog	Program Name	Phase 1	Phase 2	Phase 3	Phase 4	Phase 5	Phase 6	Duration
1	Prewash	Mains water						5'
2	Rapid 60°C	Wash in mains water at 60°C with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 60°C		55'
3	Standard 80°C	Wash in mains water at 80°C with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 75°C		62'
4	Intensive 90°C	Cold prewash in mains water	Wash in mains water at 90°C with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 75°C	65'
5	Thermal disinfection 93°C – 3 min	Thermal disinfection at 93°C for 3 min, in mains water with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 75°C		70'
6	Thermal disinfection 93°C – 10 min	Thermal disinfection at 93°C for 10 min, in mains water with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 75°C		75'
7	Intensive thermal disinfection 93°C – 10 min	Wash in mains water at 45°C	Thermal disinfection at 93°C for 10 min, in mains water with alkaline detergent	Neutralisation with acid detergent	Rinse in mains water	Rinse in demineralised water	Rinse in demineralised water at 75°C	80'
8	Service program	Charging of peristaltic pumps	Rinse in mains water					5'