



# Whatman Microbiological Monitors

Save up to 70% in  
sample preparation time.

Whatman microbiological monitors provide a complete system solution for liquid sample preparation. Each monitor is a single-use, pre-sterilized filtering unit consisting of a measured filter funnel, base, pad, membrane, removable lid and plug. With no need for flaming or sterilization, Whatman monitors cut multiple steps out of the testing process, saving your lab up to 70% in preparation time. These ready-to-use 100 mL units are suited for monitoring contaminants in all types of samples, from raw materials to finished products.

## Features and Benefits

- **Rapid testing** With no need to sterilize funnels or filter base between samples, testing time can be reduced by up to 70%.
- **No flaming required** Combined filtration unit minimizes the risk of cross-contamination.
- **All-in-one system** Filtration unit easily converts to a Petri dish, which can be labeled and incubated for culturing.
- **Reduced contamination** Single-use materials virtually eliminate cross-contamination between funnel and membrane.
- **Reproducible results** All-in-one filtration unit reduces the chance of external error.
- **Easy handling** Ready-to-use, pre-sterilized monitors are simple to use.

## Applications

Microbiological analysis of:

- Potable water
- Waste water
- Soft drinks
- Beer and wine
- Dairy products
- Pharmaceuticals and cosmetics

## Workflow



1. Sample is poured into monitor unit and filtered.



2. Funnel is removed from monitor unit.  
*(Workflow continued on rear)*

Whatman®

## Workflow (continued)



3. Microbiological media is added to membrane.



4. Lid is placed on top of membrane, forming a Petri dish.



### Whatman Quality

Whatman is a global leader in separations technology and is known in the scientific community for providing innovative Life Science products and solutions. Our instinct for simplification accelerates the rate of discovery, reduces costs and saves time. For more information, visit [www.whatman.com](http://www.whatman.com).

Whatman is a registered trademark of the Whatman Group.

**North America** Whatman Inc.  
200 Park Avenue, Suite 210  
Florham Park, NJ 07932 USA  
Tel: 1-800-WHATMAN (US and Canada)  
Fax: 1-973-245-8329  
E-mail: [info@whatman.com](mailto:info@whatman.com)

**Europe** Whatman International Ltd  
Springfield Mill, James Whatman Way  
Maidstone, Kent ME14 2LE UK  
Tel: + 44 (0)1622 676670  
Fax: + 44 (0)1622 677011  
E-mail: [information@whatman.com](mailto:information@whatman.com)

Whatman GmbH  
Hahnstrasse 3  
D-37586 Dassel  
Germany  
Tel: + 49 5561 791 0  
Fax: + 49 5561 791 536  
E-mail: [information@whatman.com](mailto:information@whatman.com)

**Japan** Whatman Japan KK  
JPR Ichigaya Building 6F  
Kudan-Minami 4-7-15, Chiyoda-ku  
Tokyo 102-0074 Japan  
Tel: +81 (0)3 5215 1242  
Fax: +81 (0)3 5215 1246  
E-mail: [japaninfo@whatman.com](mailto:japaninfo@whatman.com)

**Asia Pacific** Whatman Asia Pacific Pte Ltd  
171 Chin Swee Road  
#08-01 San Centre  
Singapore 169877  
Tel: +65 6534 0138  
Fax: +65 6534 2166  
E-mail: [wap@whatman.com](mailto:wap@whatman.com)

### WHATMAN CATALOG ORDERING INFORMATION

Product	Description	Qty/Pkg	Cat. Number	
			47 mm	56 mm
Monitor	0.2 µm, white/black grid	50	10 497 511	10 497 603
Monitor	0.45 µm, white/black grid	50	10 497 500	10 497 600
Monitor	0.45 µm, white/black grid, packaged*	50	10 497 501	—
Monitor	0.45 µm, black/white grid	50	10 497 502	10 497 601
Monitor	0.8 µm, black/white grid	50	10 497 503	10 497 602

\* Individually wrapped, sterile

# Whatman®

Leaders in Separations Technology  
[www.whatman.com](http://www.whatman.com)